MIKRO 2.0 / MIKRO 2.0 R

Inhalt des Dokuments / content of the document

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Operating instructions (EN)

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Operating instructions

MIKRO 2.0 / MIKRO 2.0 R

Translation of the original operating instructions

©2025 - All rights reserved

Andreas Hettich GmbH

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D-78532 Tuttlingen, Germany

Telephone: +49 (0)7461 705-0

Fax: +49 (0)7461 705-1125

Email: info@hettichlab.com, service@hettichlab.com

Internet: www.hettichlab.com

Table of contents

1	About this document.	6
1.1	Use of this document.	6
1.2	Gender reference.	6
1.3	Symbols and labels in this document.	6
1.4	Explanation of terms.	7
2	Safety.	7
2.1	Intended use.	7
2.2	Foreseeable misuse.	7
2.3	Use of accessories.	8
2.4	Service life.	8
2.5	Personnel qualifications.	8
2.5.1	Authorisation levels.	8
2.6	Personal protective equipment.	9
2.7	Operator obligations.	9
2.8	Safety instructions.	10
3	Device overview.	12
3.1	Technical data.	12
3.2	Rating plate.	13
3.3	Certifications and logos.	14
3.4	European registration.	15
3.5	Important labels on the packaging.	15
3.6	Important labels on the device.	16
3.7	Returns.	16
3.8	Operating and indicator elements.	17
3.8.1	Control.	17
3.8.2	Controls.	17
3.9	Scope of delivery.	18
4	Transport and storage.	18
4.1	Transport and storage conditions.	18
4.2	Transport dimensions and transport weight.	19
4.3	Fastening the transport lock.	20
5	Commissioning.	20
5.1	Unpacking the centrifuge.	20
5.2	Removing the transport lock.	20
5.3	Setting up and connecting the centrifuge.	20
5.4	Switching on for the first time.	22
5.5	Switching the centrifuge on and off.	24
6	Operation.	25
6.1	Opening and closing the lid.	25
6.2	Removing and inserting the rotor.	25
6.3	Fitting and removing the rotor lid.	26
6.4	Loading.	27

6.5	Centrifugation parameters.	29
6.5.1	Relative centrifugal force, RCF.	29
6.5.2	Centrifugation of substances or mixtures of substances with a density higher than 1.2 kg/dm ³	29
6.5.3	Ramp-up and brake levels.	30
6.6	User interface.	31
6.6.1	Dealing with messages.	31
6.6.2	Basic displays.	32
6.6.3	Keypad.	34
6.7	Menus.	35
6.7.1	Main menu.	35
6.7.2	Quick Adjust Menu.	36
6.7.3	Program menu.	37
6.7.4	Settings menu.	37
6.8	Centrifugation.	37
6.8.1	Continuous operation.	37
6.8.2	Centrifugation with time setting.	38
6.8.3	Short-term centrifugation.	39
6.9	Quick stop function.	41
6.10	Software operation.	41
6.10.1	Using Quick Adjust.	41
6.10.2	Program menu.	43
6.10.2.1	Program 'Pre-cooling'.	43
6.10.2.2	Creating a program.	44
6.10.2.3	Retrieving a program.	45
6.10.2.4	Changing a program.	45
6.10.2.5	Creating program favourites.	47
6.10.2.6	Retrieving program favourites.	48
6.10.2.7	Deleting program favourites.	49
6.10.2.8	Copying a program.	50
6.10.2.9	Renaming a program.	51
6.10.2.1	0 Deleting a program.	52
6.10.2.1	1 Starting a centrifugation run in the Program menu.	53
6.10.3	Displaying parameters of the last centrifugation run.	53
6.10.4	Settings menu.	54
6.10.4.1	Device information.	54
6.10.4.2	User settings.	57
6.10.4.3	Service settings and Advanced settings.	68
6.10.4.4	Open source licences.	68
7	Maintenance, cleaning and care.	69
7.1	Cleaning.	69
7.2	Disinfection.	70
7.3	Maintenance.	71

8	Troubleshooting.	73
8.1	Fault description.	73
8.2	Warning descriptions.	77
8.3	Performing a MAINS RESET.	81
8.4	Emergency release.	81
8.5	Replacing the mains input fuse.	82
9	Spare parts.	82
10	Disposal.	82
10.1	General instructions.	82
11	Free and Open Source Software.	84
12	Glossary.	86
13	Index.	87

1 About this document

1.1 Use of this document

- Read this document carefully and in full before commissioning the device for the first time.
Observe other enclosed instruction sheets where necessary.
- This document is part of the device and must be kept within easy reach.
- This document must be included if the device is passed on to a third party.
- The current version of the document in the available languages can be found on the manufacturer's website: ➔ <https://www.hettichlab.com/de/download-center/>

1.2 Gender reference

The employed masculine or feminine language form is to facilitate reading. In the spirit of equal treatment, corresponding terms apply in principle to all genders and do not imply any valuation.

1.3 Symbols and labels in this document

General symbols

The following markers are used in this document to highlight instructions, results, listings, references and other elements:

Marker	Explanation
1.  2.  3.  ... 	Step-by-step instructions
	Results of action steps
	References to sections of the document and other applicable documents
■ ... ■ ...	Listings without a fixed order
[Buttons]	Controls (for example: buttons, switches)
'Indicator'	Indicator elements (for example: signal lights, screen elements)

Warnings

Hazard levels	Meaning
DANGER	Indicates a hazard which, if not avoided, will result in death or serious injury.
WARNING	Indicates a hazard which, if not avoided, could result in death or serious injury.
CAUTION	Indicates a hazard which, if not avoided, could result in minor or moderate injuries.

Hazard levels	Meaning
NOTE	Indicates a hazard which, if not avoided, could result in damage to property.

Worth knowing

Recommendations	Meaning
Tip/Info	Indicates information for efficient operation and know-how for effective use of the device.

1.4 Explanation of terms

The term "centrifuge" employed in these operating instructions refers to the device referred to on the cover sheet.

2 Safety

2.1 Intended use

The centrifuge **MIKRO 2.0 General Purpose, MIKRO 2.0 R General Purpose** is a general laboratory device that is placed on the market in accordance with Machinery Directive 2006/42/EC (a changeover to Machinery Directive 2023/1230 will occur in July 2026).

The device is used to separate aqueous solutions and suspensions up to a density of 1.2 kg/dm³. The user can set the variable physical parameters within the limits specified by the device.

The centrifuge may only be used by qualified personnel in closed laboratories.

Disclaimer

Intended use also includes observing all information in the operating instructions and performance of all the required inspection and maintenance work. Any other use or use beyond this scope is considered improper. Andreas Hettich GmbH shall not be liable for any damage arising from such non-compliant use.

If the centrifuge is installed in another device or integrated into a system, the manufacturer of the system as a whole is responsible for its safety.

2.2 Foreseeable misuse

Use of the centrifuge outside the defined conditions is considered improper use.

Among other things, the following are considered improper use:

- Use of the centrifuge in explosive environments.
- Use of the centrifuge in radioactive environments.
- Use of the centrifuge in biologically contaminated environments.
- Use of the centrifuge in chemically contaminated environments.
- Use of the centrifuge in corrosive environments.
- Centrifugation of combustible materials.
- Centrifugation of explosive materials.
- Centrifugation of radioactive materials.
- Centrifugation of chemically highly reactive materials.
- Centrifugation of corrosive materials.

2.3 Use of accessories

Only use accessories approved by the manufacturer for the intended use of the centrifuge.

2.4 Service life

The planned service life of the device is 7 years.

Service life of the accessory in terms of the number of cycles and/or service life as indicated on the accessory.

2.5 Personnel qualifications

All control and safety devices must only be operated by trained persons.

Qualified mechanics

For the purposes of this documentation, qualified mechanics are persons who are familiar with the design, mechanical installation, commissioning, troubleshooting and maintenance of the product and who have the following qualifications:

- Qualification / training in the field of mechanics in accordance with national regulations.

Qualified electricians

For the purposes of this documentation, qualified electricians are persons who are familiar with the design, electrical installation, commissioning, troubleshooting and maintenance of the product and who have the following qualifications:

- Qualification / training in electrical engineering in accordance with national regulations.

Qualified refrigeration technicians

For the purposes of this documentation, qualified refrigeration technicians are persons who are familiar with the design and installation of refrigeration and chilling equipment, and the commissioning, troubleshooting and maintenance of the product and who have the following qualifications:

- Qualification / training in the field of refrigeration and cooling technology in accordance with national regulations.
- An appropriate certificate of competence (e.g. in accordance with ISO 22712) for working with A3 refrigerants must be available.

Trained person

For the purposes of this documentation, these are persons who have received instruction enabling them to carry out work in the areas of transportation, storage and operation.

2.5.1 Authorisation levels

Authorisation level 1

"Restricted rights"

A user who uses the system to complete a task. Centrifugations can be started and stopped. Changing programs or viewing and changing system settings is not permitted.

If you select authorisation level 1, you can specify which programs can be executed.

The user can only use some functions after a user with authorisation level 3 has entered a PIN.

Example: loader / assistant / student

Authorisation level 2

"Standard rights"

Authorisation level 1 rights are included.

A user with advanced knowledge. The user is allowed to change the value settings in Quick Adjust. This has a corresponding effect on short-term centrifugation. Event messages/logs can be viewed. Readout and analysis of system information is possible to a limited extent.

Some functions can only be used by authorisation level 2 users after an authorisation level 3 user has entered a PIN.

Authorisation level 3

"Administrator"

Authorisation level 2 rights are included.

A user who is allowed to make changes to the system and programs. Readout and analysis of system information is possible.

The user must set the PIN when using the device for the first time.

Example: laboratory manager / product manager

2.6 Personal protective equipment

Observe national and local regulations on safety and accident prevention. Select and provide protective clothing such as protective footwear, safety gloves and protective goggles in alignment with the respective activities and the expected risks.

2.7 Operator obligations

Ensure that all the work listed below is carried out only by qualified personnel:

- Transportation, unpacking, lifting
- Setup
- Operation
- Maintenance
- Disposal

The operator must ensure that persons working on the device observe the rules and regulations as well as the following instructions:

- National and local regulations pertaining to safety, accident prevention and environmental protection.
- Do not install or commission any damaged devices.
- The necessary protective equipment must be provided.
- Operate the device only when it is in perfect working order.
- Periodically check that personnel are working safely and in compliance with the safety information in the operating instructions.
- Check protective equipment regularly to ensure it is serviceable.
- Do not remove attached safety and warning notices and ensure they remain legible.
- Persons under the influence of alcohol, drugs or medicines that affect their ability to react must not operate or maintain the device.
- Keep the operating instructions available in the place where the device is used.

- Pass the operating instructions on to the buyer when the device is sold.

2.8 Safety instructions

Reporting of events and incidents

In the event of incidents or notifiable events involving the device or its accessories, these must be reported to the manufacturer and, where applicable, to the competent authority where the user and/or the patient is registered.

Manufacturer:

Andreas Hettich GmbH

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78532 Tuttlingen, Germany

Phone: +49 7461 705 0

E-mail: info@hettichlab.com

DANGER

Risk of contamination

Inadequate cleaning or failure to observe the cleaning instructions can lead to contamination risks.

- Observe national and local regulations on safety and accident prevention.
- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Comply with laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

DANGER

Risk of fire and explosion

Risk of accidents, injuries or damage to property due to fire or explosion.

- Observe regulations and directives for handling chemicals and hazardous substances.
- Do not use corrosive chemicals.
- Do not use dangerous chemicals.
- Do not use corrosive extraction agents.
- Do not use strong acids.

WARNING

Risk of injury

Inadequate or late maintenance can result in injuries.

- Observe maintenance intervals.
- Check the device for visible damage or defects.
If any visible damage or defects are present, immediately remove the device from service and inform a service technician.

**! WARNING****Risk of electric shock**

Liquids that penetrate the device may cause electric shocks.

- The device must be protected from external contact with liquids.
- Do not pour any liquids inside the device.
- The original transport packaging must be used when transporting the device.

**! WARNING****Contamination with hazardous substances and substance mixtures**

Observe the following actions for substances and substance mixtures that are toxic, radioactive and/or contaminated with pathogenic microorganisms:

- As a rule, use only centrifuge tubes with special screw caps for hazardous substances.
- Use sealable centrifuge tubes with a biosafety system for materials of risk groups 3 and 4.
- If no biosafety system is used, the device is not micro-biologically tight in the sense of standard EN / IEC 61010-2-020.
- Contact the manufacturer if necessary.

**WARNING****Risk of injury and device damage due to the rotor**

A loose rotor can cause injuries and device damage.

- Check that the rotor is firmly seated.
- Follow maintenance intervals.

**CAUTION****Risk of injury**

Long hair and clothing can get caught in the rotor during manual movement.

- Tie long hair back.
- Do not allow garments to hang in the centrifuging chamber.

**NOTICE****Damage**

Incorrect voltage or frequency.

- Only operate the device according to the specifications on the rating plate.
- Compliance with the instructions for use.

**NOTICE****Damage**

Aborting the program early can cause damage to the device and samples.

- Do not switch off, perform an emergency release or pull out the mains plug.

3 Device overview

3.1 Technical data

Manufacturer	Andreas Hettich GmbH D-78532 Tuttlingen	
Model	MIKRO 2.0	MIKRO 2.0 R
Type	2440	2460
Mains voltage (±10%)	100-240 V 1 ~	100-240 V 1 ~
Mains frequency	50-60 Hz	50-60 Hz
power consumption	270 VA	420 VA
Refrigerant	-	R290
Refrigerant content	-	35 g
Global warming potential (GWP)	-	0,02
max. capacity	12 x 5 ml	
max. permissible density	1.2 kg/dm³	
max. speed (RPM)	16100	
max. acceleration (RCF)	25000	
max. kinetic energy	6400 Nm	
Obligation to perform checks (DGUV Rules 100-500) (valid only in Germany)	No	
Ambient conditions (EN / IEC 61010-1):		
Installation site	indoors only	
Altitude	up to 3000 m above sea level	
Ambient temperature	2 °C to 40 °C	5 °C to 35 °C
Humidity	maximum relative humidity 80% for temperatures up to 31 °C, decreasing linearly to 50% relative humidity at 40 °C.	

Overvoltage category (IEC 60364-4-443)	II	
Pollution level	2	
Device protection class	I not suitable for use in potentially explosive atmospheres.	
EMC:		
Emitted EM interference, EM interference immunity	EN / IEC 61326-1 Class B FCC Class B	EN / IEC 61326-1 Class B FCC Class B
Noise level (rotor-dependent)	≤58 dB(A)	≤61 dB(A)
Dimensions:		
Width	240 mm	240 mm
Depth	389 mm	538 mm
Altitude	225 mm	250 mm
Weight	approx. 18 kg	approx. 28 kg

3.2 Rating plate

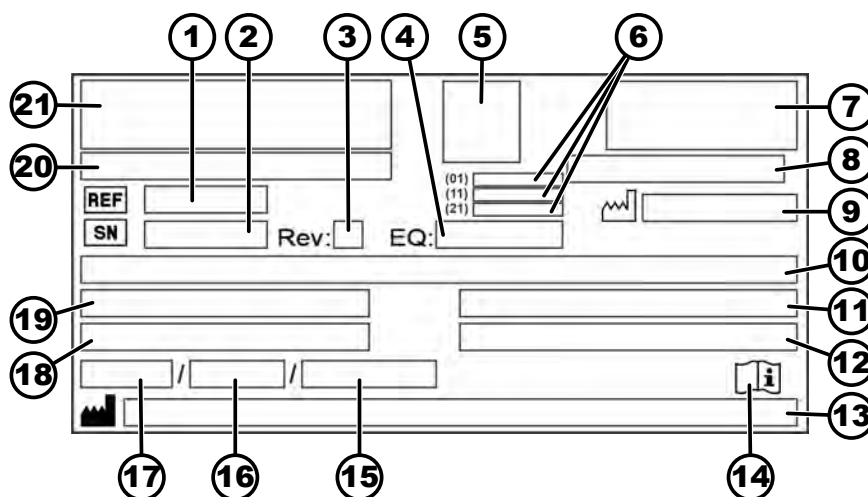
Fig. 1: Rating plate

1	REF Item number	12	Maximum permissible density
2	SN Serial number	13	Manufacturer's address
3	Revision	14	Follow the user manual
4	Equipment number	15	Maximum permissible pressure (PS) in the refrigeration system (only for devices with cooling)


5	UDI code	16	Refrigerant content (only for devices with cooling)
6	Content of the UDI code: (01) Global Trade Item Number (GTIN) (11) Date of manufacture (21) Serial number	17	Refrigerant type (only for devices with cooling)
7	EAC mark, CE mark	18	Maximum speed in revs per minute
8	Country of manufacture	19	Current draw (A) or power consumption (VA)
9	 Date of manufacture	20	Device designation
10	Mains voltage and mains frequency	21	Manufacturer's logo
11	Maximum kinetic energy		

3.3 Certifications and logos

Certifications

 ISO 9001 Certified Quality Management System www.tuvsud.com/ms-cert	ISO 9001 Quality management system in accordance with ISO 9001
 ISO 14001 Certified Environmental Management System www.tuvsud.com/ms-cert	ISO 14001 Environmental management in accordance with ISO 14001
 Certified Quality System ISO 13485 mdc	EN ISO 13485 Quality management in accordance with ISO 13485

Logos

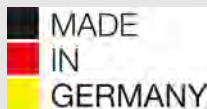 MADE IN GERMANY	Made in Germany Device was developed and made in Germany.
---	---


3.4 European registration

Device conformity

Device conformity according to EU directives.

3.5 Important labels on the packaging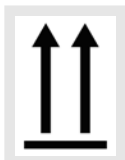

TOP

This is the correct upright position of the shipping container for transport and/or storage.


FRAGILE GOODS

The contents of the shipping container are fragile, so it must be handled with care.


PROTECT FROM MOISTURE

Protect the transport packaging from moisture and keep it in a dry environment.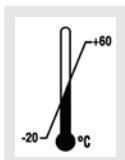

TEMPERATURE LIMITATION

The shipping container must be stored, transported and handled within the indicated temperature range (-20 °C to +60 °C).

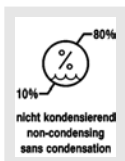
HUMIDITY LIMITATION

The shipping container must be stored, transported and handled within the indicated air humidity range (10 % to 80 %, non-condensing).

STACK LIMITATION BASED ON QUANTITY

Maximum number of identical packages that may be stacked on the lowest package, "n" standing for the number of packages allowed. The lowest package is not included in "n".

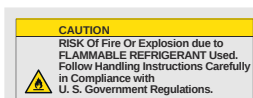Warning of risk of fire and explosion due to flammable refrigerant

3.6 Important labels on the device

The signs on the device must not be removed or covered, or have anything pasted over them.

Attention, general danger area.


Ensure you read the instructions for commissioning and operation and observe the safety instructions before using the device.

Biohazard warning.

Warning of fire hazard.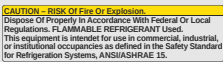

This warning indicates the risk of fire or explosion caused by refrigerants.

Direction of rotation of the rotor.

The orientation of the arrow indicates the rotor's direction of rotation.

Direction of rotation of the emergency release.

Symbol for the separate collection of electrical and electronic equipment, in accordance with Directive 2012/19/EU (WEEE).

Use in European Union countries, Norway and Switzerland.

3.7 Returns

An original Return Material Authorisation (RMA) form from the manufacturer must always be requested for a return. Secure and reliable acceptance and booking in of the goods with the manufacturer is not possible without an original RMA form from the manufacturer. The Return Material Authorisation (RMA) form contains a Declaration of No Objection (UBE), which must be completed in full and enclosed with the return.

If a leak is suspected or confirmed by the service department, the refrigeration circuit must be evacuated prior to return. The same applies in the event of a crash because it is not certain that all the refrigerant will escape.

If the device and/or accessories are returned to the manufacturer, the complete return shipment must be cleaned and decontaminated by the sender. If returns are not cleaned and/or decontaminated or are insufficiently cleaned and/or decontaminated, this will be performed by the manufacturer and charged to the sender.

The original transport locks must be attached for return shipment, see ➔ Chapter 4 'Transport and storage' on page 18. The device must be shipped in its original packaging.

3.8 Operating and indicator elements

3.8.1 Control

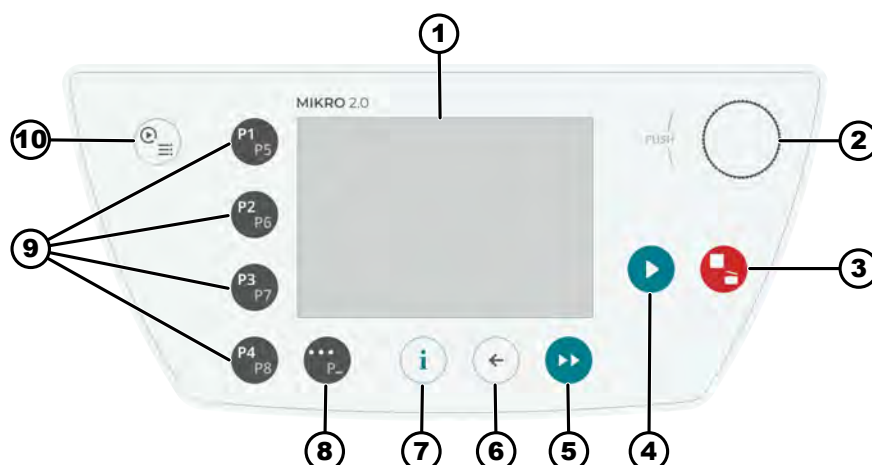
Fig. 2: Control (device without cooling)

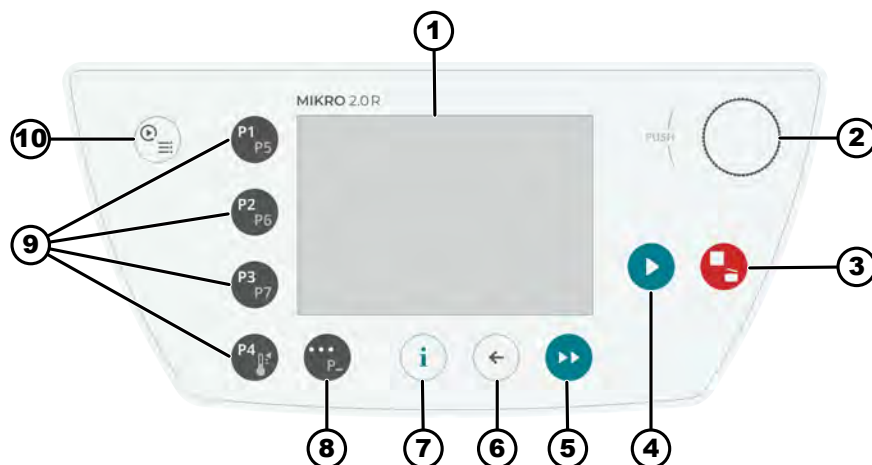
Fig. 3: Control (device with cooling)

3.8.2 Controls

Number	Button and display elements	Description
1	Screen	Displays, information and menus
2	Rotary, push button	- Rotate: set the parameters and change to the next menu item. - Press: jump to the next parameter and select a menu item.

Number	Button and display elements	Description
3	STOP, OPEN	- Press once: centrifuge brakes with set braking ramp. - Press twice, or press and hold (quick stop): centrifuge brakes as quickly as possible. - Press when the centrifuge is at a standstill: open the lid.
4	START	Start centrifugation run. A centrifugation run can be started from Quick Adjust or from the program.
5	SHORT-TERM CENTRIFUGATION	The centrifugation run takes place as long as the button is being pressed.
6	BACK	- Back to the menus. - Save changed values.
7	INFO	Call up operational data and rotor data.
8	CONTEXT BUTTON, TOGGLE BUTTON	- You can access the context menu by pressing the button as soon as '...' is displayed in a menu. - To select the second level of favourite programs, hold down the button and simultaneously press the desired program shortcut button.
9	SHORTCUT BUTTONS	- Shortcut buttons for favourite programs P1-P8. - Shortcut buttons for favourite programs P1-P7 for MIKRO 2.0 R only. - Temperature control program for MIKRO 2.0 R only.
10	QUICK ADJUST, PROGRAM	- Quick Adjust: Quick setting of parameters. The values can subsequently be transferred to a program. - Program: Call up and edit programs.

3.9 Scope of delivery

- 1 offset screwdriver
- 1 power cable
- 1 user manual
- For products distributed in Switzerland:
1 Import Switzerland sticker

4 Transport and storage

4.1 Transport and storage conditions

Transport conditions

NOTICE

The centrifuge's refrigeration circuit contains flammable refrigerant.

- Comply with possible transport restrictions (e.g. according to ADR or IATA DGR).

NOTICE

Damage

The device may be damaged if it is not secured during transport.

- Secure the transport locks before transport.
- Observe the transport instructions.

NOTICE

Danger due to condensation in the event of temperature differences

Moisture may damage electrical components.

- Ensure that all surfaces are dry prior to commissioning or maintenance.
- If the temperature changes, wait until the device or component has become acclimatised.
- Prevent moisture from penetrating into sensitive components.
- If moisture forms, switch off the device immediately and allow it to dry properly.

- Before transporting, fasten the transport lock and disconnect the device from the mains socket.
- The transport temperature must be between -20 °C and +60 °C.
- Humidity must not be condensing. Humidity must be between 10% and 80%.
- Be aware of the weight of the device.
- When transporting using a transport aid (e.g., a pallet truck), the transport aid must be able to carry at least 1.6 times the transport weight of the device.
- Secure the device to prevent it tipping over and falling down during transport.
- Never transport the device sideways or upside down.

Storage conditions

- The device must be stored in the original packaging.
- Only store the device in dry rooms.
- The same restriction regarding room size is applicable to storage as well as operation. ➔ *Chapter 2.8 'Safety instructions' on page 10.*
- The storage temperature must be between -20 °C and +60 °C.
- Humidity must not be condensing. Humidity must be between 10% and 80%.

4.2 Transport dimensions and transport weight

Article	Packaging dimensions in mm (L x W x H)	Weight in kg
MIKRO 2.0 / MIKRO 2.0 R	240 x 389 x 255 / 240 x 538 x 250	approx. 20 kg / approx. 33 kg

4.3 Fastening the transport lock

Personnel:

- Trained user

1. ➤ Open the lid.
2. ➤ Insert the foam cylinder.
3. ➤ Close the lid.

5 Commissioning

5.1 Unpacking the centrifuge

CAUTION

Danger of crushing due to parts falling out of the transport packaging.

- Keep the device balanced during the unpacking process.
- Only open the packaging at the points provided for this purpose.

CAUTION

Risk of injury from lifting heavy loads.

- Provide an adequate number of helpers.
- Note the weight. See ➔ *Chapter 3.1 'Technical data' on page 12.*

NOTICE

Damage to the device due to improper lifting.

- Do not lift the centrifuge by the control panel or the control panel holder.

5.2 Removing the transport lock

Personnel:

- Trained user

1. ➤ Open the lid.
2. ➤ Remove the foam cylinder.
3. ➤ Store the foam cylinder in a safe place.
4. ➤ Close the lid.

5.3 Setting up and connecting the centrifuge

Setting up the centrifuge

WARNING

Operation of the unit in an explosive, corrosive or radioactively contaminated atmosphere is not in accordance with the intended use.

**WARNING****Risk of injury**

Due to insufficient clearance from the centrifuge.

- As per EN / IEC 61010-2-020, no persons, hazardous materials or objects may be present within a safety zone of 300 mm around the centrifuge during a centrifugation run.
- Maintain a distance of 300 mm from the ventilation slots and ventilation openings of the centrifuge.
- The centrifuge's ventilation openings must never be blocked.

**CAUTION****Risk of crushing and damage**

Changes in position caused by vibration may cause the device to fall.

- Place the device on a stable and level surface.
- Select an installation area that is capable of bearing the weight of the device.
- Observe national and local regulations on safety and accident prevention.

**NOTICE****Damage**

Any deviation from the external temperature conditions will result in damage to the samples and device.

- Comply with the maximum and minimum permissible ambient temperatures.
- Avoid placing the device next to heat sources.
- Do not place the device in direct sunlight.
- Protect the device from frost.
- Maintain the required clearance around the device.

Personnel:

- Trained user

1. ➤ Place the device on a stable and level surface.
2. ➤ Maintain a distance of 300 mm around the device.
3. ➤ Comply with the ambient conditions in the technical data (➔ *Chapter 3.1 'Technical data' on page 12*).

Connecting the centrifuge

**NOTICE****Damage**

Property damage caused by unauthorised personnel.

- Do not allow persons without the appropriate authorisation to perform any work on or make changes to devices.
- Only authorised personnel may carry out maintenance and repairs.
- Obtain the manufacturer's approval or guidance before carrying out any work on the device.

**NOTICE****Danger due to condensation in the event of temperature differences**

Moisture may damage electrical components.

- Ensure that all surfaces are dry prior to commissioning or maintenance.
- If the temperature changes, wait until the device or component has become acclimatised.
- Prevent moisture from penetrating into sensitive components.
- If moisture forms, switch off the device immediately and allow it to dry properly.

Personnel:

- Trained user

1. ➤ A type B residual current circuit breaker must be used if the device is additionally protected with a residual current circuit breaker in the building installation.

When using a different type, the residual current circuit breaker may either not switch off the unit if there is a fault on the unit, or it may switch off the unit even though there is no fault on the unit.

2. ➤ Check whether the mains voltage and mains frequency match the specification on the rating plate.
3. ➤ Connect the device to a standard mains socket using the mains cable.

5.4 Switching on for the first time

Personnel:

- Authorisation level 3

1. ➤ Set the mains switch to */I/*.

➡ Progress bar is displayed.

The company logo is briefly displayed.

Animation and introductory text are briefly displayed.

'Initial start-up' is displayed.

'Start configuration' is highlighted in blue.

2. ➤ Press *[Rotary, push button]*.

➡ Language selection is displayed.

3. ➤ Rotate *[Rotary, push button]* until the desired language is highlighted in blue.
4. ➤ Press *[Rotary, push button]*.
 - The date format and date selection are displayed.
5. ➤ Rotate *[Rotary, push button]* until the desired date format setting is selected.
6. ➤ Press *[Rotary, push button]*.
 - The day setting is highlighted in blue and displayed in bold.
7. ➤ Rotate *[Rotary, push button]* until the desired value for the day is selected.
8. ➤ Press *[Rotary, push button]*.
 - The month setting is highlighted in blue and displayed in bold.
9. ➤ Rotate *[Rotary, push button]* until the desired value for the month is selected.
10. ➤ Press *[Rotary, push button]*.
 - The year setting is highlighted in blue and displayed in bold.
11. ➤ Rotate *[Rotary, push button]* until the desired value for the year is selected.
12. ➤ Press *[Rotary, push button]*.
 - The time format and time selection are displayed.
13. ➤ Rotate *[Rotary, push button]* until the desired time format setting is selected.
14. ➤ Press *[Rotary, push button]*.
 - The hour setting is highlighted in blue and displayed in bold.
15. ➤ Rotate *[Rotary, push button]* until the desired value for the hours is selected.
16. ➤ Press *[Rotary, push button]*.
 - The minute setting is highlighted in blue and displayed in bold.
17. ➤ Rotate *[Rotary, push button]* until the desired value for the minutes is selected.
18. ➤ Press *[Rotary, push button]*.
 - The PIN input menu is displayed.
19. ➤ Rotate *[Rotary, push button]* to select the desired number, then press *[Rotary, push button]* to enter the number.

To delete a number, rotate *[Rotary, push button]* until the  symbol is highlighted in blue, then press *[Rotary, push button]*.

Repeat the process until all 8 numbers have been entered.

- ✓ symbol is highlighted in blue.
20. ➤ Press *[Rotary, push button]*.
 - The PIN confirmation menu is displayed.

21. ➤ Rotate [Rotary, push button] to select the desired number, then press [Rotary, push button] to enter the number.

To delete a number, rotate [Rotary, push button] until the  symbol is highlighted in blue, then press [Rotary, push button].

Repeat the process until all 8 numbers have been entered.

- ✓ symbol is highlighted in blue.

22. ➤ Press [Rotary, push button].

- The settings are stored.

'Commissioning successful' is displayed.

'Restart centrifuge' is highlighted in blue.

23. ➤ Press [Rotary, push button].

- Centrifuge is restarted.

5.5 Switching the centrifuge on and off.

Switching the centrifuge on

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

- Set the mains switch to [I].

- The following displays appear one after the other:

- Progress bar
- Company logo

- Quick Adjust menu is displayed.

The last menu used before the device was switched off is always displayed. If a program was the last thing to be executed, for example, the Program menu is also displayed when the device is switched on.

To simplify and clarify presentation, in this document it is always assumed that the Quick Adjust menu is started.

Switching off the centrifuge

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is stationary.

- Set the mains switch to [O].

- Device shuts down.

Blue background with symbol is displayed.

6 Operation

6.1 Opening and closing the lid

Opening the lid

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The centrifuge is switched on.

The rotor is stationary.

→ Press the *[STOP, OPEN]* button.

- ➡ The lid unlocks by means of a motor.

Closing the lid

CAUTION

Crushing hazard when closing the lid.

Danger of fingers getting crushed when the closing motor pulls the lid against the seal.

- No parts of the operator's body should be in the hazard zone of the lid when closing the lid.
- To close the lid, press on the lid from above.

NOTICE

Damage to the device caused by the lid slamming.

- Close the lid slowly.
- Do not slam the lid.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

→ Close the lid and press the front edge of the lid down gently.

- ➡ The lid locks using a motor.

6.2 Removing and inserting the rotor

Rotor removal

Personnel:

- Trained user

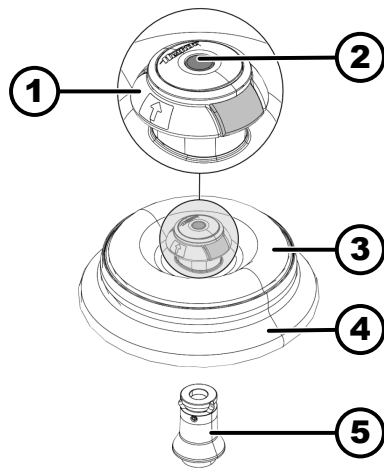
Fig. 4: Inserting and removing the rotor

- 1 Removal aid
- 2 Release pin
- 3 Rotor lid
- 4 Rotor
- 5 Rotor mount

1. ➤ Open the centrifuge lid.
2. ➤ Grip beneath the removal aid (1) with your index finger and middle finger.
3. ➤ Press the release pin (2) down with your thumb.
4. ➤ Pull the removal aid (1) upwards.
5. ➤ Lift the rotor (4) with the rotor lid (3) upwards from the rotor mount (5) and remove.

Inserting the rotor

Personnel:

- Trained user

The centrifuge lid is open.

1. ➤ Clean the rotor mount (5) and rotor hole.
2. ➤ Grip beneath the removal aid (1) with your index finger and middle finger.
 - ➡ Removal aid (1) moves upwards.
3. ➤ Place the rotor (4) vertically on the rotor mount (5).
4. ➤ Check whether the rotor (4) is correctly inserted.

6.3 Fitting and removing the rotor lid

Removing the rotor lid

Personnel:

- Trained user

The centrifuge lid is open.

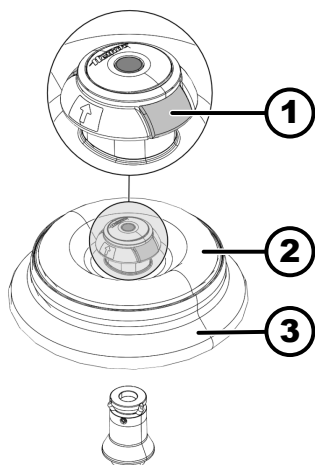
Fig. 5: Fitting and removing the rotor lid

- 1 Lid release
- 2 Rotor lid
- 3 Rotor

Fitting the rotor lid

Personnel:

- Trained user

The centrifuge lid is open.

1. Press the lid release (1) together with your index finger and thumb.
2. Place the rotor lid (2) vertically on the rotor (3).
3. Check whether the rotor lid (2) is correctly engaged.
Check whether the lid release (1) is in the correct position.

6.4 Loading

Filling centrifuge tubes

WARNING

Risk of injury from contaminated sample material

Leakage of contaminated sample material from the sample tube during centrifugation.

- When working with hazardous substances, only use centrifuge tubes with special screw caps.
- For materials in risk groups 3 and 4, additionally use a biosafety system in accordance with the WHO 'Laboratory Biosafety Manual' alongside sealable centrifuge tubes.

NOTICE

Damage

Device damage due to corrosive substances.

- Do not centrifuge highly corrosive substances.
- Do not allow corrosive substances to enter the device.
- Check the device regularly for corrosion.

Standard glass centrifuge tubes can be loaded up to RCF 4000 (DIN 58970 part 2).

Personnel:

- Trained user

→ Fill centrifuge tubes outside the centrifuge.

The maximum capacity of the centrifuge tubes specified by the manufacturer must not be exceeded.

With angle rotors, the centrifuge tubes must only be filled to the extent that no liquid can be ejected from the tubes during the centrifugation run.

It must be ensured that there is a uniform fill level in the tubes in order to keep the weight differences in the centrifuge tubes as low as possible.

Symbols on the angle rotor

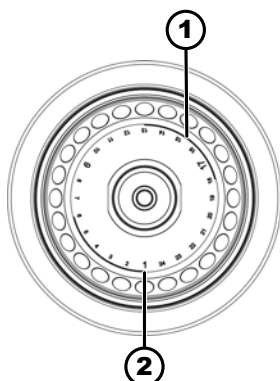
Fig. 6: Rotor with symbols

- 1 Division of the rotor into four segments
- 2 Division of the rotor into three segments


The following symbols appear on the rotor to help with even rotor loading:

- Division of the rotor into four segments.
Two thick arcs are used for marking.
Loading aid for even distribution of the centrifugation tubes over four rotor segments.
- Division of the rotor into three segments.
Location numbers are used for marking, and these appear in a larger font.
Loading aid for even distribution of the centrifugation tubes over three rotor segments.

Loading the angle rotors

Personnel:

- Trained user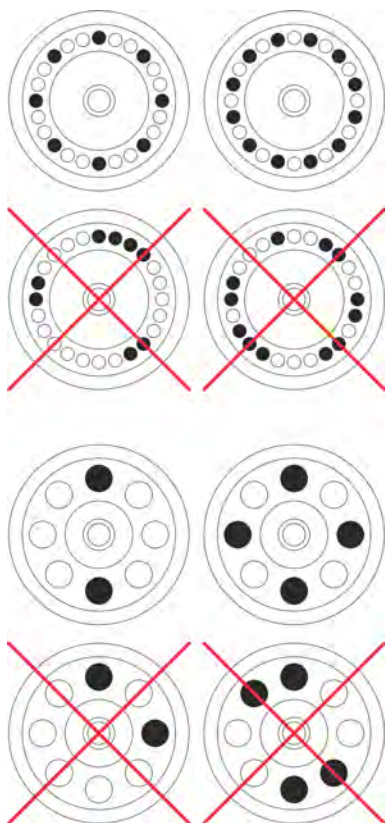

1. ➤ Check that the rotor is firmly seated.
2. ➤ The centrifuge tubes must be distributed evenly over all locations on the rotor.

No liquid must be allowed to enter the rotor and the centrifuging chamber when loading the rotor.

With rotors, the centrifuge tubes must only be filled to the extent that no liquid can be ejected from the tubes during the centrifugation run.

The weight of the permissible filling capacity is indicated on each rotor. The weight must not be exceeded.

6.5 Centrifugation parameters

6.5.1 Relative centrifugal force, RCF

The relative centrifugal force (RCF) is dependent on the speed and the centrifuging radius.

RCF is specified as a multiple of the acceleration due to gravity (g).

The RCF is a dimensionless numerical value and is used to compare the separation and sedimentation performance.

$$RCF = \left(\frac{RPM}{1000} \right)^2 * r * 1,118$$

$$RPM = \sqrt{\frac{RCF}{r * 1,118}} * 1000$$

RCF = Relative Centrifugal Force

RPM = revolutions per minute, $[RPM] = \text{min}^{-1}$.

r = distance from the axis of rotation to the bottom of the centrifuge tube, $[r] = \text{mm}$.

6.5.2 Centrifugation of substances or mixtures of substances with a density higher than 1.2 kg/dm³

The density of the substances or mixtures of substances must not exceed 1.2 kg/dm³ during centrifugation at maximum speed. The speed must be reduced for substances or substance mixtures with a higher density. The permissible speed can be calculated using the following formula:

$$\text{Reduced speed } (n_{red}) = \sqrt{\frac{1,2}{\text{Greater density } [\text{kg/dm}^3]}} * \text{maximum speed } [RPM]$$

For example: Maximum speed 4000 RPM, density 1.6 kg/dm³

$$n_{red} = \sqrt{\frac{1,2(\text{kg/dm}^3)}{1,6(\text{kg/dm}^3)}} * 4000 \text{ RPM} = 3464 \text{ RPM}$$

If, in exceptional cases, the maximum load indicated on the bucket is exceeded, the speed must also be reduced. The permissible speed can be calculated using the following formula:

$$\text{Reduced speed } (n_{red}) = \sqrt{\frac{\text{maximum load [g]}}{\text{actual load [g]}}} * \text{maximum speed [RPM]}$$

For example: Maximum speed 4000 RPM, maximum load 300 g, actual load 350 g

$$n_{red} = \sqrt{\frac{300 \text{ g}}{350 \text{ g}}} * 4000 \text{ RPM} = 3703 \text{ RPM}$$

Please contact the manufacturer if you are not sure.

6.5.3 Ramp-up and brake levels

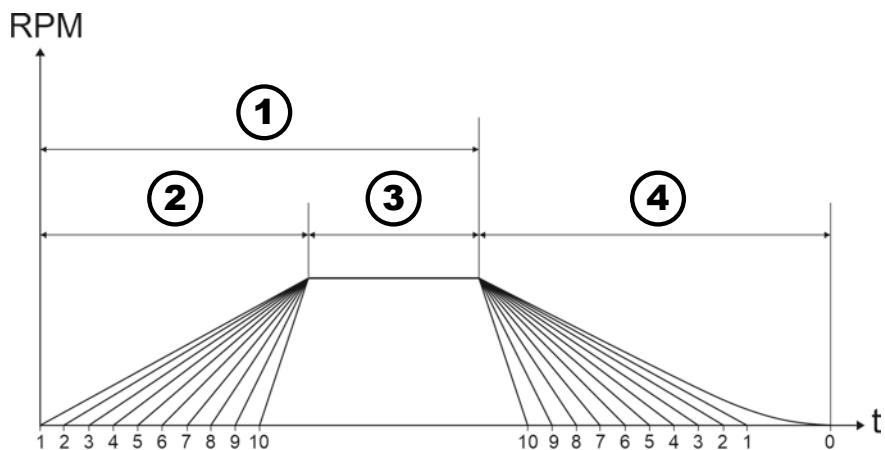
Fig. 7: Ramp-up and brake levels

- 1 Runtime (runtime measurement after the start of the centrifugation run)
- 2 Ramp-up levels
- 3 Runtime (runtime measurement after reaching the speed)
- 4 Brake levels

Runtime

You can set how the runtime is measured:

- Runtime measurement starts after the start of the centrifugation run (1).
- Runtime measurement starts after reaching centrifugation speed (3).

To set the runtime measurement, see ➔ Chapter 6.10.1 'Using Quick Adjust' on page 41.

Ramp-up levels

10 ramp-up levels can be set (2).

- Ramp-up level 1: longest time until the speed is reached.
- Ramp-up level 10: shortest time until the speed is reached.

To set the ramp-up levels, see ➔ Chapter 6.10.1 'Using Quick Adjust' on page 41.

Brake levels

11 brake levels can be set (4).

- Brake level 10: shortest time until the rotor comes to a standstill.
- Brake level 1: second longest time until the rotor comes to a standstill.
- Brake level 0: unbraked ramp-down of the rotor. Longest time until the rotor comes to a standstill.

To set the brake levels, see ➔ *Chapter 6.10.1 'Using Quick Adjust' on page 41.*

6.6 User interface

6.6.1 Dealing with messages

An information, warning or error message is displayed in the information line of the user interface if the centrifuge is in a state that the user needs to be informed about.

For certain messages, both an audible signal (see ➔ *Chapter 6.10.4.2.9 'Audible signal' on page 60*) and a visual signal (see ➔ *Chapter 6.10.4.2.8 'Visual signal' on page 60*) can also be set.

Warning messages can only be acknowledged when a user with authorisation level 3 has entered a PIN.

Information about the centrifugation run

- Show details

Information about the centrifugation run appears in the information line.


1. ➤ Press the *[INFO]* button.
 - ➔ Details of the information about the centrifugation run are displayed on the user interface.
2. ➤ Press the *[BACK]* button.
 - ➔ Details of the information about the centrifugation run are hidden.

- Acknowledge information

Information about the centrifugation run appears in the information line.

- Press *[Rotary, push button]* or the *[BACK]* button, or open the lid.
 - ➔ Information is acknowledged.

Warning messages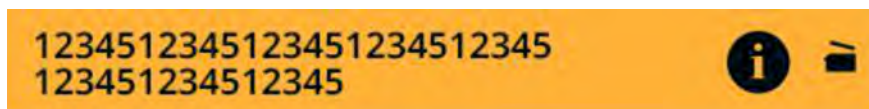

- Show details

Warning message appears in the information line.

1. ➤ Press the *[INFO]* button.
 - ➔ Details of the warning message are displayed on the user interface.
2. ➤ Press the *[BACK]* button.
 - ➔ Details of the warning message are hidden.

- Acknowledge message


Warning message appears in the information line.

1. ➤ *[Rotary, push button]* or press the *[BACK]* button.
 - ➔ Query is displayed.
2. ➤ Press *[Rotary, push button]*.
 - ➔ Message is acknowledged.

A PIN must be entered for authorisation level 1.

It may be necessary to enter a PIN for authorisation level 2.

Error messages

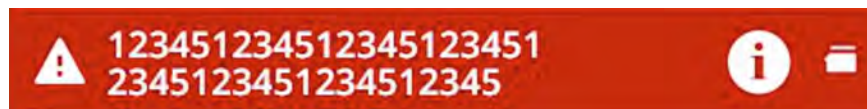- Show details

Error message appears in the information line.

1. ➤ Press the **[INFO]** button.
 - Details of the error message are displayed on the user interface.
2. ➤ Press the **[BACK]** button.
 - Details of the error message are hidden.

- Acknowledge message

Error message appears in the information line.

- Switching the centrifuge off and on again, see ➔ *Chapter 5.5 'Switching the centrifuge on and off.'* on page 24.
- If the error message persists, contact the manufacturer's service department.

6.6.2 Basic displays

All illustrated user interfaces and their contents are for illustration purposes only.

The parameters of the actual programs may differ from those shown here.

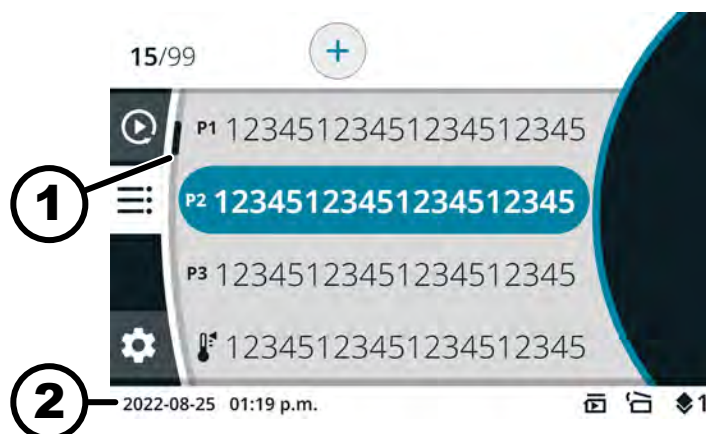
Fig. 8: Program menu

- 1 Scroll bar
- 2 Status bar

Scroll bar

In the various menus, the scroll bars indicate the length of the list and the position of the selected entry.

Status bar

The following symbols can be displayed in the status bar:

Symbol	Meaning
	Indicates the progress of the centrifugation run.
	Start of the centrifugation run as soon as the lid is closed.

Symbol	Meaning
	For uncooled centrifuges only: Lid opens automatically after a successful centrifugation run
	Authorisation level Level 1: Restricted rights Level 2: Standard rights Level 3: Administrator
	The set parameters do not match the rotor used.

Information lines

Fig. 9: Information line

1 Information line

The following symbols can be displayed in the information lines:

Symbol	Meaning
	The lid is closed.
	The lid is open.
	More information can be displayed by pressing the [INFO] button.
	error message
	Wait
	Short-term centrifugation

Symbol	Meaning
	Copy
	Settings
	Action performed successfully.
	Action cancelled.

6.6.3 Keypad

Fig. 10: Keypad

- 1 Country-specific keypad
- 2 Delete input
- 3 Space bar
- 4 Toggle between upper and lower case letters
- 5 Special characters
- 6 Letters and numbers

Letters and numbers

- 1. ➤ Rotate [Rotary, push button] until the desired letter or number is highlighted in blue.
- 2. ➤ Press [Rotary, push button].
 - The letter or number is transferred to the input field.

Space bar

- 1. ➤ Rotate [Rotary, push button] until the space bar (3) is highlighted in blue.
- 2. ➤ Press [Rotary, push button].
 - A space is transferred to the input field.

Toggle between upper and lower case letters

- 1. ➤ Rotate [Rotary, push button] until the symbol (4) is highlighted in blue.
- 2. ➤ Press [Rotary, push button].
 - Upper or lower case letters are displayed.

Special characters

1. ➤ Rotate *[Rotary, push button]* until the symbol (5) is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - The special characters are displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired special character is highlighted in blue.
4. ➤ Press *[Rotary, push button]*.
 - The special character is transferred to the input field.

Delete input

1. ➤ Rotate *[Rotary, push button]* until the symbol (2) is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - The last entry is deleted from the input field.

Country-specific keypad

1. ➤ Rotate *[Rotary, push button]* until the language code of the country (1) is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - The country-specific keypad for the displayed country is displayed.
 - The next possible language code (1) is displayed in alphabetical order.

6.7 Menus

6.7.1 Main menu

Fig. 11: Main menu

- 1 Quick Adjust menu
- 2 Program menu
- 3 Settings menu
- 4 Status bar

Selecting a menu

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

1. ➤ Rotate *[Rotary, push button]* until the desired menu symbol is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - The background of the selected menu symbol turns white.
 - Menu contents are displayed.

Changing the menu

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

A menu is selected.

1. ➤ Press the *[BACK]* button.
 - The background of the menu symbol is highlighted in blue.
 - The main menu is displayed.
2. ➤ Rotate *[Rotary, push button]* until the desired menu symbol is highlighted in blue.
3. ➤ Press *[Rotary, push button]*.
 - The background of the selected menu symbol turns white.
 - Menu contents are displayed.

Procedure if you only want to toggle between the Quick Adjust menu and the Program menu:

- Press the *[QUICK ADJUST, PROGRAM]* button.
 - You can toggle back and forth between the Quick Adjust menu and the Program menu.
 - Menu contents are displayed.

6.7.2 Quick Adjust Menu

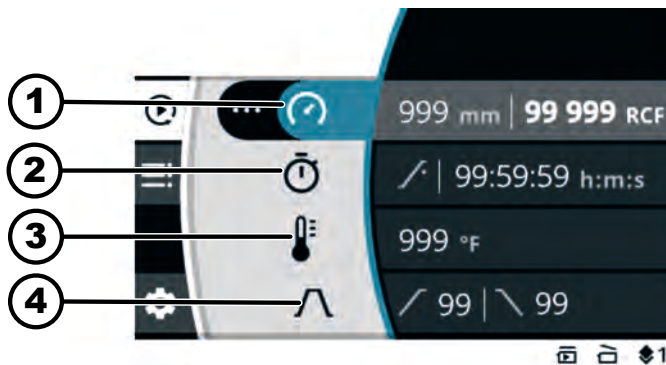
Fig. 12: Quick AdjustMenu

- 1 Centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 2 Runtime
- 3 Temperature (only for devices with cooling)
- 4 Ramp-up and brake level

To select the Quick Adjust menu, see ➔ *Chapter 6.7.1 'Main menu' on page 35*

To set the parameters, see ➔ *Chapter 6.10.1 'Using Quick Adjust' on page 41*

6.7.3 Program menu

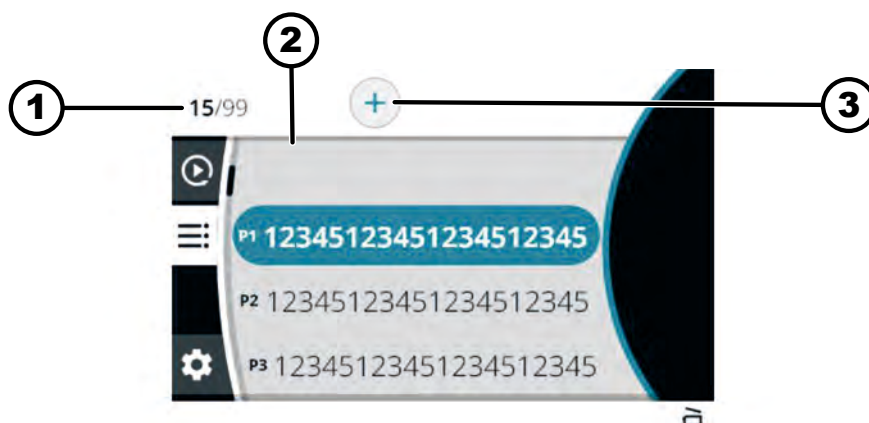
Fig. 13: Program menu

- 1 Number of programs created / maximum number of possible programmes
- 2 List of existing programs
- 3 Create new program

Menu selection, see ➔ Chapter 6.7.1 'Main menu' on page 35

For further settings, see ➔ Chapter 6.10.2 'Program menu' on page 43

6.7.4 Settings menu

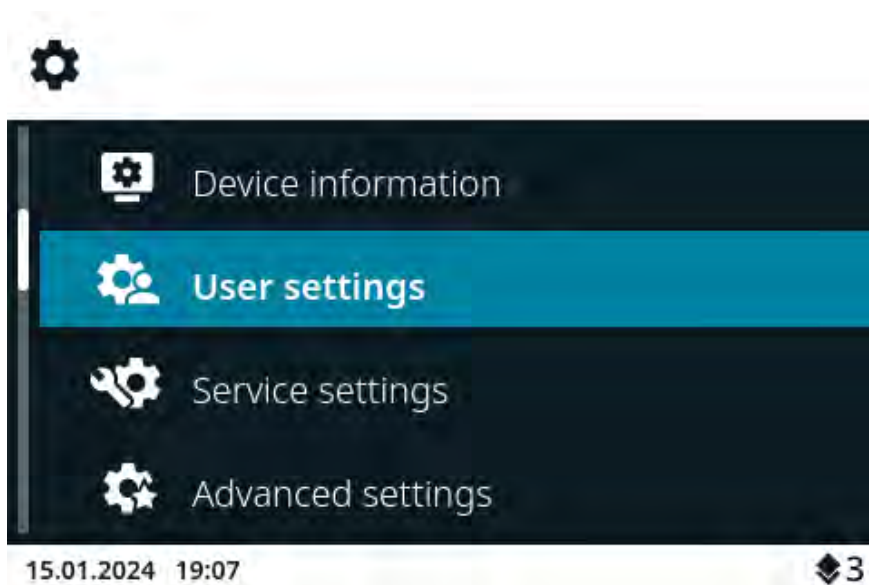
Fig. 14: Settings menu

To select the Settings menu, see ➔ Chapter 6.7.1 'Main menu' on page 35

For further settings, see ➔ Chapter 6.10.4 'Settings menu' on page 54

6.8 Centrifugation

6.8.1 Continuous operation

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3


The rotor is installed and the centrifuge lid is closed.

The  symbol is set as the runtime in the Quick Adjust menu or the Program menu of the selected program. ➔ *Chapter 6.10.1 'Using Quick Adjust' on page 41*

1. ➔ Press the **[START]** button.

➔ The centrifugation run is started.

The following summary is displayed during the centrifugation run: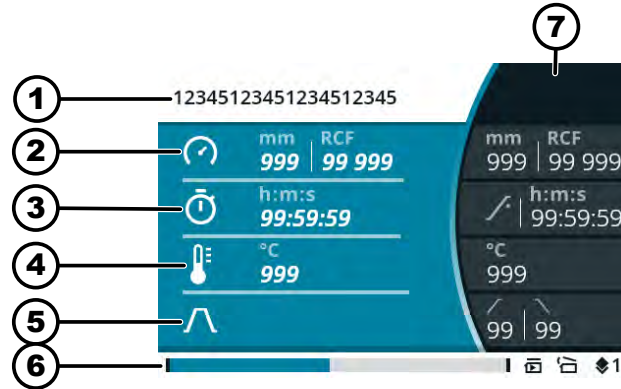

- 1 Selected program or Quick Adjust
- 2 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 3 Current runtime
- 4 Current temperature (for devices with cooling)
- 5 Symbol for ramp-up and brake level
- 6 Status bar
- 7 Set target values

2. ➔ Press the **[STOP, OPEN]** button to end the centrifugation run.

➔ Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

3. ➔ If required: Press the **[INFO]** button after the centrifugation run.

➔ Centrifugation run parameters are displayed.

4. ➔ If required: Press the **[BACK]** button.

➔ Centrifugation run parameters are hidden.

5. ➔ Opening the lid

or

Press the **[BACK]** button.

➔ The original screen is displayed.

6.8.2 Centrifugation with time setting

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3


The rotor is installed and the centrifuge lid is closed.

The desired parameters are set in the Quick Adjust menu or the program selected from the Program menu. ➔ *Chapter 6.10.1 'Using Quick Adjust' on page 41* ➔ *Chapter 6.7.3 'Program menu' on page 37*

1. ➤ Press the *[START]* button.

➔ The centrifugation run is started.

The following summary is displayed during the centrifugation run: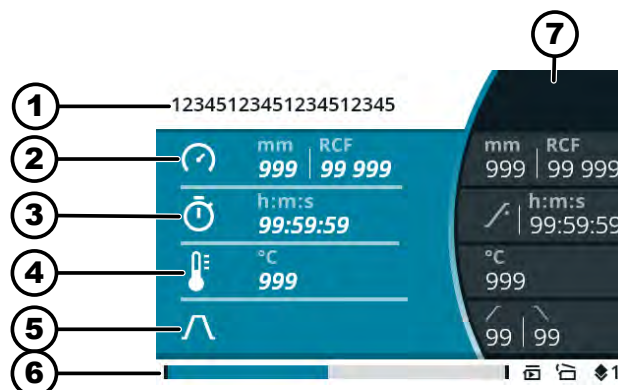

- 1 Selected program or Quick Adjust
- 2 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 3 Remaining runtime
- 4 Current temperature (for devices with cooling)
- 5 Symbol for ramp-up and brake level
- 6 Status bar
- 7 Set target values

2. ➤ Wait until the runtime has elapsed, or press the *[STOP, OPEN]* button to interrupt the centrifugation run.

➔ Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

3. ➤ If required: Press the *[INFO]* button after the centrifugation run.

➔ Centrifugation run parameters are displayed.

4. ➤ If required: Press the *[BACK]* button.

➔ Centrifugation run parameters are hidden.

5. ➤ Opening the lid

or

Press the *[BACK]* button.

➔ The original screen is displayed.

6.8.3 Short-term centrifugation

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

The desired parameters are set in the Quick Adjust menu. → Chapter 6.10.1 'Using Quick Adjust' on page 41


1. ➤ Press and hold the [SHORT-TERM CENTRIFUGATION] button.

➡ The centrifugation run is started.

Short-term centrifugation can be started when the Quick Adjust menu, the main menu or the program list is displayed.

The following summary is displayed during the centrifugation run: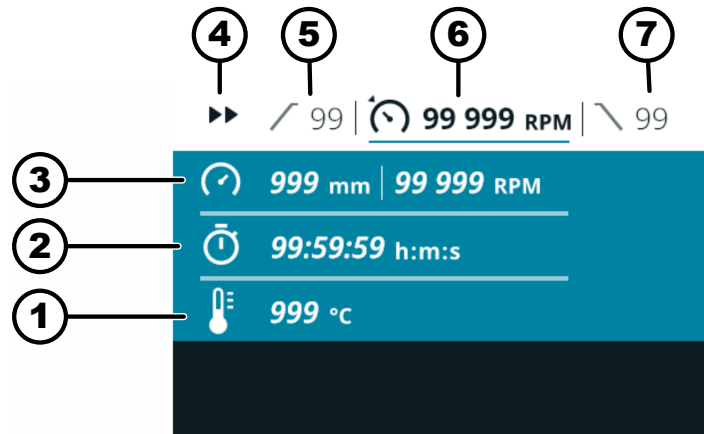

- 1 Current temperature (for devices with cooling)
- 2 Current runtime
- 3 Current centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 4 Short-term centrifugation indicator
- 5 Set ramp-up level
- 6 Set speed (RPM) or relative centrifugal force (RCF)
- 7 Set brake level

2. ➤ Release the [SHORT-TERM CENTRIFUGATION] button to end the centrifugation run.

➡ Ramp-down takes place with the set brake level.

The lid can be opened shortly after the rotor has come to a standstill.

Information line with information about the centrifugation run is displayed.

3. ➤ If required: Press the [INFO] button.

➡ Centrifugation run parameters are displayed.

4. ➤ If required: Press the [BACK] button.

➡ Centrifugation run parameters are hidden.

5. ➤ Opening the lid

or

Press the [BACK] button.

➡ The original screen is displayed.

6.9 Quick stop function

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

➞ Press the *[STOP, OPEN]* button twice
or

Press and hold the *[STOP, OPEN]* button.

➞ Ramp-down takes place with the fastest possible braking.

6.10 Software operation

6.10.1 Using Quick Adjust

Quick Adjust menu is selected, ➞ Chapter 6.7.2 'Quick Adjust Menu' on page 36.

Personnel:

- Authorisation level 2
- Authorisation level 3

Change parameters while at a standstill

The following description assumes that the selection 'RPM/RCF' is set as 'Starting point in Quick Adjust' in the 'User settings' menu. Otherwise, setting of the runtime commences. ➞ Chapter 6.10.4.2.1 'Starting point in Quick Adjust' on page 57

Changed values are applied directly.

1. ➞ Rotate *[Rotary, push button]* to change the speed (RPM) or relative centrifugal force (RCF).

2. ➞ Press *[Rotary, push button]*.

➞ Value is saved.

The next menu symbol is highlighted in blue.

The runtime is displayed in bold and can be changed.

3. ➞ Rotate *[Rotary, push button]* to change the runtime.

For centrifuging in continuous operation, set the  symbol.

The symbol is located between '99:59:59 h:m:s' and '00:00:01 h:m:s'.

4. ➞ Press *[Rotary, push button]*.

➞ Value is saved.

The next menu symbol is highlighted in blue.

For centrifuges with cooling: The temperature is displayed in bold and can be changed.

For centrifuges without cooling: The ramp-up level is displayed in bold and can be changed. The next two steps are skipped.

5. ➞ Rotate *[Rotary, push button]* to change the temperature.

Changing the relative centrifugal force (RCF), centrifuging radius

6. ➤ Press *[Rotary, push button]*.
 - Value is saved.
 - The symbol for the ramp-up and brake level is highlighted in blue.
 - The ramp-up level is displayed in bold and can be changed.
7. ➤ Rotate *[Rotary, push button]* to change the ramp-up level.
8. ➤ Press *[Rotary, push button]*.
 - The brake level is displayed in bold and can be changed.
9. ➤ Rotate *[Rotary, push button]* to change the brake level.

Personnel:

- Authorisation level 2
- Authorisation level 3

Menu symbol for the speed (RPM), relative centrifugal force (RCF) and the centrifuging radius is selected.

1. ➤ Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
2. ➤ Rotate *[Rotary, push button]* to select 'RPM' or 'RCF'.
 - Selected value is displayed in bold.
3. ➤ Press *[Rotary, push button]*.
 - Radius is selected and displayed in bold.
4. ➤ Rotate *[Rotary, push button]* to change the radius.
5. ➤ Press the *[BACK]* button.
 - The changes are stored.

Changing the runtime starting time

Personnel:

- Authorisation level 2
- Authorisation level 3

Menu symbol for the runtime is selected.

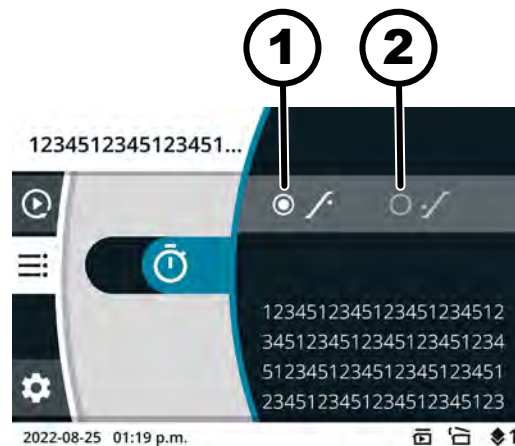
Fig. 15: Starting time - Runtime

- 1 Timing of the runtime starts after the set speed is reached
 - 2 Runtime begins after the start of the centrifugation run
1. ➤ Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
 2. ➤ Rotate *[Rotary, push button]* to change the value.
 - Selected value is displayed in bold.
 3. ➤ Press the *[BACK]* button.
 - The changes are stored.

Start centrifugation run

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

➞ Press the *[START]* button.

Changing parameters during the centrifugation run

Personnel:

- Authorisation level 2
- Authorisation level 3

Centrifugation run in the Quick Adjust menu is active.

The runtime display is set to 'classic'.

1. ➞ Press and hold the *[Rotary, push button]* button for 5 seconds.
 - ➞ The speed (RPM) or relative centrifugal force (RCF) parameters of the current centrifugation run is shown in bold in the Quick Adjust menu.
2. ➞ Rotate *[Rotary, push button]* to change the value.
3. ➞ Press *[Rotary, push button]*.
 - ➞ The runtime parameter of the current centrifugation run is shown in bold in the Quick Adjust menu.
4. ➞ Rotate *[Rotary, push button]* to change the value.
 - ➞ The next two steps are skipped for centrifuges without cooling.
5. ➞ Press *[Rotary, push button]*.
 - ➞ The temperature parameter of the current centrifugation run is shown in bold in the Quick Adjust menu.
6. ➞ Rotate *[Rotary, push button]* to change the value.
7. ➞ Press *[START]* to save the changes to the parameters in Quick Adjust or
Press *[BACK]* to discard the changes to the parameters in Quick Adjust.

6.10.2 Program menu

6.10.2.1 Program 'Pre-cooling'

(only for devices with cooling)

The 'Pre-cooling' program is used to cool the unloaded rotor to the desired temperature for the subsequent centrifugation.

Only the temperature can be set. All other parameters are fixed and cannot be changed.

The set temperature must be lower than the current temperature (actual temperature), otherwise the centrifugation run cannot be started.

The centrifugation run for pre-cooling the rotor stops automatically as soon as the set temperature is reached.

The 'Pre-cooling' program cannot be renamed or deleted.

Retrieving the program 'Pre-cooling' using the shortcut button

Personnel:

- Authorisation level 2
- Authorisation level 3

Main menu or Quick Adjust menu or Program menu is selected, ➔ *Chapter 6.7 'Menus' on page 35*

The Pre-cooling program is located in favourite place 1.

1. ➔ Press and hold [CONTEXT BUTTON, TOGGLE BUTTON].
2. ➔ Press [SHORTCUT BUTTON] of favourite place 1.
 - ➔ Program 'Pre-cooling' is highlighted in blue.

Program 'Pre-cooling' parameters are displayed.

Program menu is selected. ➔ *Chapter 6.7.3 'Program menu' on page 37*

Retrieving program 'Pre-cooling' with [Rotary, push button]

1. ➔ Rotate [Rotary, push button] until program 'Pre-cooling' is highlighted in blue.
2. ➔ Press [Rotary, push button].
 - ➔ Program parameters are displayed.

Personnel:

- Authorisation level 3

Program 'Pre-cooling' is retrieved. ➔ *'Retrieving the program 'Pre-cooling' using the shortcut button' on page 44.* ➔ *'Retrieving program 'Pre-cooling' with [Rotary, push button]' on page 44*

Changing program 'Pre-cooling'

1. ➔ Press [Rotary, push button].
 - ➔ The temperature is displayed in bold and can be changed.
2. ➔ Rotate [Rotary, push button] to change the temperature.
3. ➔ Press the [BACK] button.
 - ➔ The change is stored.

6.10.2.2 Creating a program

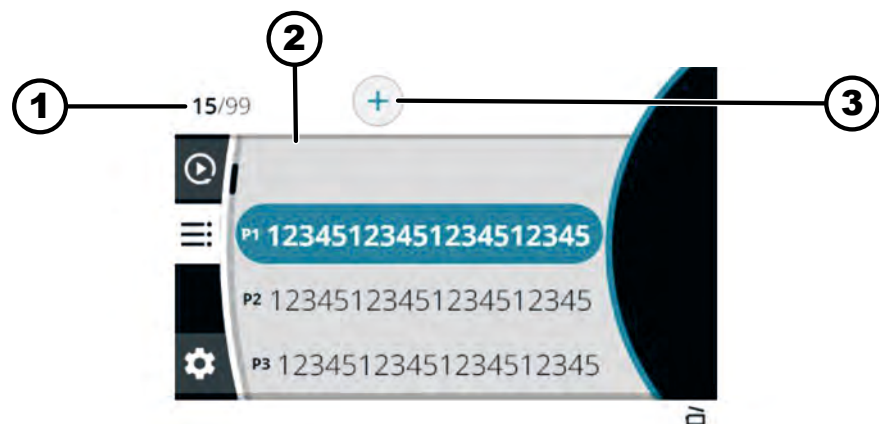
Fig. 16: Program menu

- 1 Number of programs created / maximum number of possible programmes
- 2 List of existing programs
- 3 Create new program

Personnel:

- Authorisation level 3

Program menu is selected. ➔ *Chapter 6.7.3 'Program menu' on page 37*

1. ➤ Rotate *[Rotary, push button]* to the right until the create new program symbol is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Keyboard opens.
3. ➤ Rotate and press *[Rotary, push button]* to enter the program name.

To delete a number, rotate *[Rotary, push button]* until the symbol is highlighted in blue, then press *[Rotary, push button]*.

Repeat the process until the program name has been entered.

4. ➤ Press the *[BACK]* button.
 - ✓ symbol is highlighted in blue.
5. ➤ Press *[Rotary, push button]*.
 - Program is saved with the parameters from the Quick Adjust menu.

6.10.2.3 Retrieving a program

Personnel:

- Authorisation level 2
- Authorisation level 3

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37

1. ➤ Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Program parameters are displayed.

It is still possible to select a program by rotating the *[Rotary, push button]*.
3. ➤ Press *[Rotary, push button]*.
 - Program parameters are retrieved.

6.10.2.4 Changing a program

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37
Program is called.

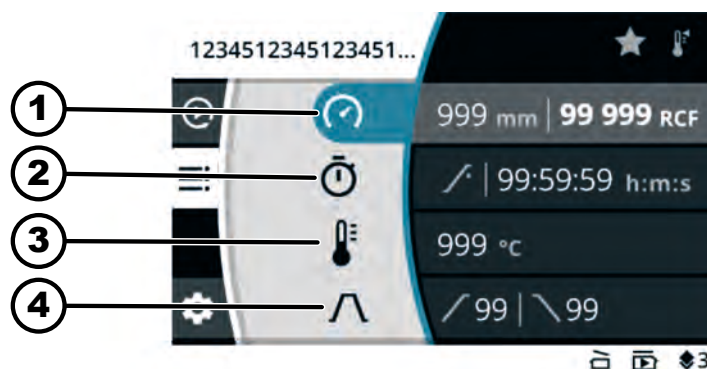
Fig. 17: Program mode menu - Change program

- 1 Centrifuging radius, speed (RPM) or relative centrifugal force (RCF)
- 2 Runtime

- 3 Temperature (only for devices with cooling)
- 4 Ramp-up and brake level

Personnel:

- Authorisation level 2
- Authorisation level 3

1.  Rotate *[Rotary, push button]* to change the speed (RPM) or relative centrifugal force (RCF).
2.  Press *[Rotary, push button]*.
 - ➡ Value is saved.
 - The next menu symbol is highlighted in blue.
 - The runtime is displayed in bold and can be changed.
3.  Rotate *[Rotary, push button]* to change the runtime.
 - For centrifuging in continuous operation, set the  symbol.
 - The symbol is located between '99:59:59 h:m:s' and '00:00:01 h:m:s'.
4.  Press *[Rotary, push button]*.
 - ➡ Value is saved.
 - The next menu symbol is highlighted in blue.
 - For centrifuges with cooling: The temperature is displayed in bold and can be changed.
 - For centrifuges without cooling: The ramp-up level is displayed in bold and can be changed. The next two steps are skipped.
5.  Rotate *[Rotary, push button]* to change the temperature.
6.  Press *[Rotary, push button]*.
 - ➡ Value is saved.
 - The symbol for the ramp-up and brake level is highlighted in blue.
 - The ramp-up level is displayed in bold and can be changed.
7.  Rotate *[Rotary, push button]* to change the ramp-up level.
8.  Press *[Rotary, push button]*.
 - ➡ The brake level is displayed in bold and can be changed.
9.  Rotate *[Rotary, push button]* to change the brake level.
10.  Press the *[BACK]* button.
 - ➡ The changes are stored.

Changing the relative centrifugal force (RCF), centrifuging radius

Personnel:

- Authorisation level 2
- Authorisation level 3

Menu symbol for the speed (RPM), relative centrifugal force (RCF) and the centrifuging radius is selected.

1.  Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
2.  Rotate *[Rotary, push button]* to select 'RPM' or 'RCF'.
 - ➡ Selected value is displayed in bold.
3.  Press *[Rotary, push button]*.
 - ➡ Radius is selected and displayed in bold.
4.  Rotate *[Rotary, push button]* to change the radius.

Changing the runtime starting time

5. Press the *[BACK]* button.
 - ➡ The changes are stored.

Personnel:

- Authorisation level 2
- Authorisation level 3

Menu symbol for the runtime is selected.

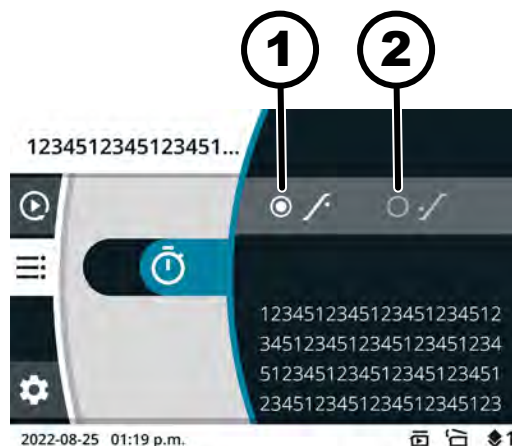
Fig. 18: Starting time - Runtime

- 1 Timing of the runtime starts after the set speed is reached
 - 2 Runtime begins after the start of the centrifugation run
1. ➤ Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
 2. ➤ Rotate *[Rotary, push button]* to change the value.
 - ➔ Selected value is displayed in bold.
 3. ➤ Press the *[BACK]* button.
 - ➔ The changes are stored.

6.10.2.5 Creating program favourites

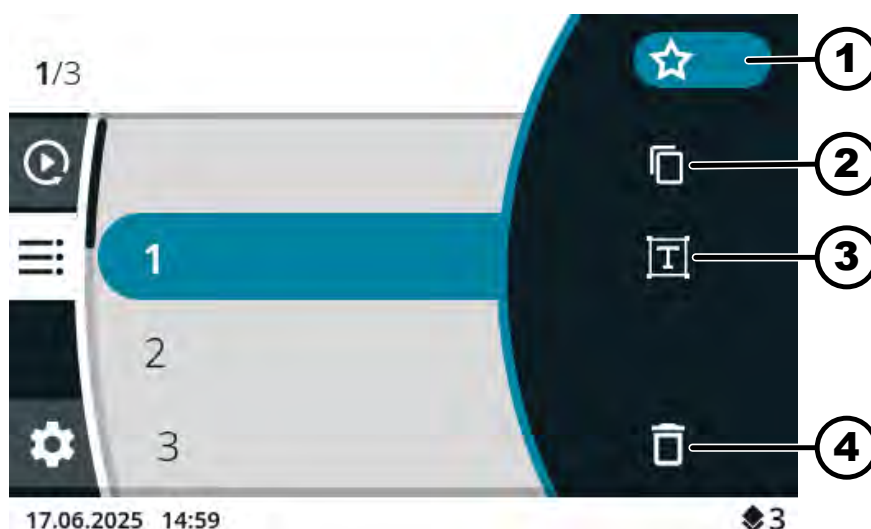
Fig. 19: Program mode symbols

- 1 "Favourites" symbol
- 2 "Copy program" symbol
- 3 "Change program name" symbol
- 4 "Delete program" symbol

Personnel:

- Authorisation level 3

Program menu is selected. ➔ *Chapter 6.7.3 'Program menu' on page 37*

1. ➤ Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Program parameters are displayed.
3. ➤ Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
 - Submenu appears.
4. ➤ Rotate *[Rotary, push button]* until the "Favourites" symbol is highlighted in blue.
5. ➤ Press *[Rotary, push button]*.
 - The Favourites submenu is displayed.
 - Favourite places that are already assigned are marked with a solid star.
6. ➤ Rotate *[Rotary, push button]* until the desired free favourite place is highlighted in blue.
7. ➤ Press *[Rotary, push button]*.
 - Program is assigned to the favourite place.

6.10.2.6 Retrieving program favourites

Personnel:

- Authorisation level 2
- Authorisation level 3

Main menu or Quick Adjust menu or Program menu is selected, ➔ *Chapter 6.7 'Menus' on page 35*

Favourite places P1 to P4

- Press the *[SHORTCUT BUTTON]* of the desired favourite place.
 - Favourite program is highlighted in blue.
 - Program parameters are displayed.

Favourite places P5 to P8

1. ➤ Press and hold *[CONTEXT BUTTON, TOGGLE BUTTON]*.
2. ➤ Press the *[SHORTCUT BUTTON]* of the desired favourite place.
 - Favourite program is highlighted in blue.
 - Program parameters are displayed.

6.10.2.7 Deleting program favourites

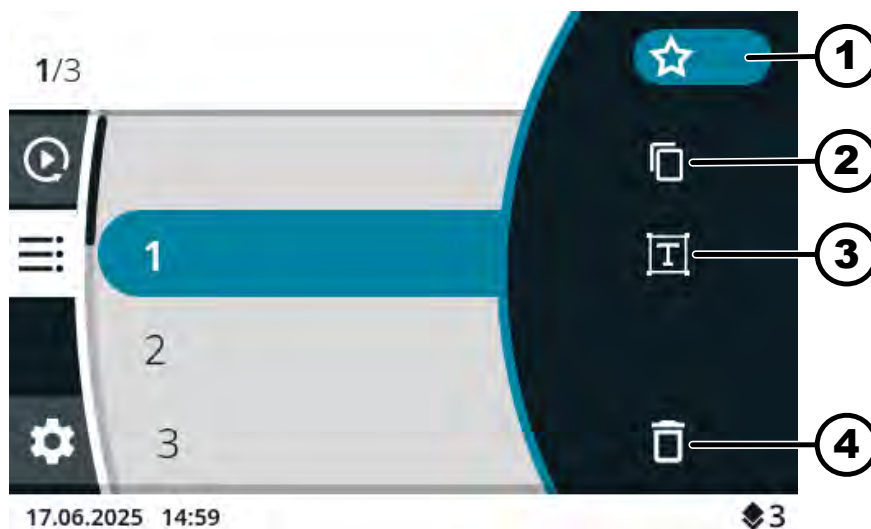
Fig. 20: Program mode symbols

- 1 "Favourites" symbol
- 2 "Copy program" symbol
- 3 "Change program name" symbol
- 4 "Delete program" symbol

Personnel:

- Authorisation level 3

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37

1. ➤ Rotate [Rotary, push button] until the desired favourite program is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Program parameters are displayed.
3. ➤ Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
 - Submenu appears.
4. ➤ Rotate [Rotary, push button] until the "Favourites" symbol is highlighted in blue.
5. ➤ Press [Rotary, push button].
 - The Favourites submenu is displayed.
 - The assigned favourite place is highlighted in blue.
6. ➤ Press [Rotary, push button].
 - Program is removed from the favourite place.

6.10.2.8 Copying a program

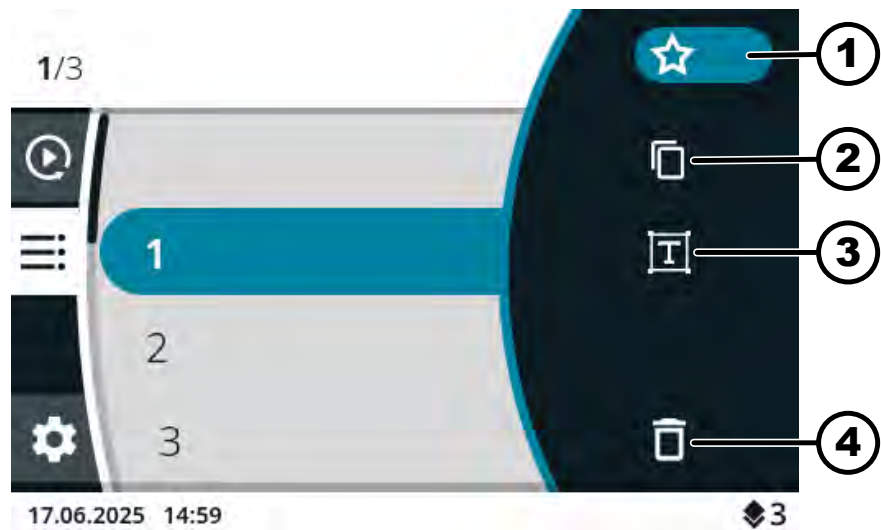
Fig. 21: Program mode symbols

- 1 "Favourites" symbol
- 2 "Copy program" symbol
- 3 "Change program name" symbol
- 4 "Delete program" symbol

Personnel:

- Authorisation level 3

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37

1. ➔ Rotate [Rotary, push button] until the desired program is highlighted in blue.
2. ➔ Press [Rotary, push button].
 - ➔ Program parameters are displayed.
3. ➔ Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
 - ➔ Submenu appears.
4. ➔ Rotate [Rotary, push button] until the "Copy program" symbol is highlighted in blue.
5. ➔ Press [Rotary, push button].
 - ➔ Keyboard opens.
6. ➔ Rotate [Rotary, push button] until the  symbol is highlighted in blue.
7. ➔ Press [Rotary, push button] repeatedly until the program name is deleted.
8. ➔ Rotate and press [Rotary, push button] to enter the program name.

To delete a number, rotate [Rotary, push button] until the  symbol is highlighted in blue, then press [Rotary, push button].

Repeat the process until the program name has been entered.

9. ➔ Press the [BACK] button.
 - ➔ ✓ symbol is highlighted in blue.

10. ➤ Press [Rotary, push button].
 - Program is saved with the copied parameters.
 - Information line indicates successful execution of the action.

6.10.2.9 Renaming a program

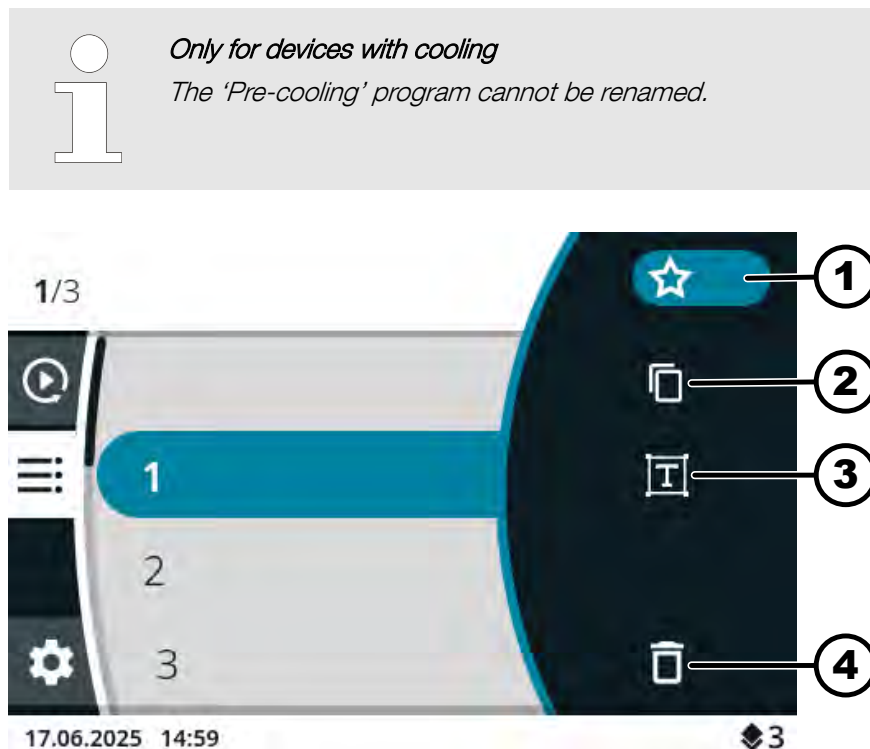
Fig. 22: Program mode symbols

- 1 "Favourites" symbol
- 2 "Copy program" symbol
- 3 "Change program name" symbol
- 4 "Delete program" symbol

Personnel:

- Authorisation level 3

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37

1. ➤ Rotate [Rotary, push button] until the desired program is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Program parameters are displayed.
3. ➤ Press the [CONTEXT BUTTON, TOGGLE BUTTON] button.
 - Submenu appears.
4. ➤ Rotate [Rotary, push button] until the "Change program name" symbol is highlighted in blue.
5. ➤ Press [Rotary, push button].
 - Keyboard opens.
6. ➤ Rotate [Rotary, push button] until the  symbol is highlighted in blue.
7. ➤ Press [Rotary, push button] repeatedly until the program name is deleted.

8. ➤ Rotate and press *[Rotary, push button]* to enter the program name.

*To delete a number, rotate *[Rotary, push button]* until the  symbol is highlighted in blue, then press *[Rotary, push button]*.*

Repeat the process until the program name has been entered.

9. ➤ Press the *[BACK]* button.
- ✓ symbol is highlighted in blue.
10. ➤ Press *[Rotary, push button]*.
- Program is saved with the new name.

6.10.2.10 Deleting a program

Only for devices with cooling

The program 'Pre-cooling' cannot be deleted.

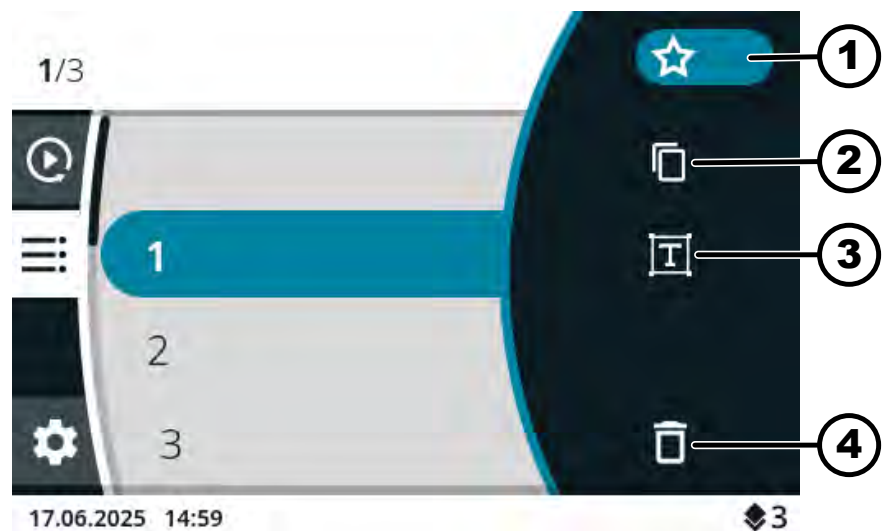
Fig. 23: Program mode symbols

- 1 "Favourites" symbol
- 2 "Copy program" symbol
- 3 "Change program name" symbol
- 4 "Delete program" symbol

Personnel:

- Authorisation level 3

Program menu is selected. ➔ Chapter 6.7.3 'Program menu' on page 37

1. ➤ Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
- Program parameters are displayed.
3. ➤ Press the *[CONTEXT BUTTON, TOGGLE BUTTON]* button.
- Submenu appears.

4. ➤ Rotate *[Rotary, push button]* until the "Delete program" symbol is highlighted in blue.
5. ➤ Press *[Rotary, push button]*.
 - Confirmation query appears
6. ➤ Rotate *[Rotary, push button]* until the ✓ symbol is highlighted in blue.
7. ➤ Press *[Rotary, push button]*.
 - Program is deleted.

6.10.2.11 Starting a centrifugation run in the Program menu

Starting a centrifugation run from the program list

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

Program menu is selected. The program names are displayed. ➔ *Chapter 6.7.3 'Program menu' on page 37*

1. ➤ Rotate *[Rotary, push button]* until the desired program is highlighted in blue.
2. ➤ Press the *[START]* button.

Starting a centrifugation run using favourite program

Personnel:

- Authorisation level 2
- Authorisation level 3

The rotor is installed and the centrifuge lid is closed.

Main menu or Quick Adjust menu or Program menu is selected, ➔ *Chapter 6.7 'Menus' on page 35*

1. ➤ Favourite places P1 to P4

Press the *[SHORTCUT BUTTON]* of the desired favourite place.

Favourite places P5 to P8

Press and hold *[CONTEXT BUTTON, TOGGLE BUTTON]*.

Press the *[SHORTCUT BUTTON]* of the desired favourite place.

- The program names are displayed.
- Favourite program is highlighted in blue.

2. ➤ Press the *[START]* button.

6.10.3 Displaying parameters of the last centrifugation run

Displaying parameters of the last centrifugation run

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

1. ➤ Press the *[INFO]* button.
 - The parameters of the last centrifugation run are displayed in the Quick Adjust menu, the program list (only when the program names are displayed) and the main menu.
2. ➤ Press the *[BACK]* button or the *[Rotary, push button]*.
 - Parameter display is ended.

6.10.4 Settings menu

Selection of submenus

1. ➤ Rotate *[Rotary, push button]* until the desired menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.

For further settings, see:

➤ Chapter 6.10.4.1 'Device information' on page 54

➤ Chapter 6.10.4.2 'User settings' on page 57

➤ Chapter 6.10.4.3 'Service settings and Advanced settings' on page 68

➤ Chapter 6.10.4.4 'Open source licences' on page 68

6.10.4.1 Device information

6.10.4.1.1 System information

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'System information' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.

The following information can be found in the menu:

- Device name
- Sales number
- Equipment number
- Serial number
- HBI software version
- CPU serial number
- ZS software version
- Manufacturer URL
- Next service
- Date/time UTC

6.10.4.1.2 Rotor information

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Rotor info' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired rotor is highlighted in blue.

4. ➤ Press *[Rotary, push button]*.

➤ Menu is displayed.

The following information can be found in the menu:

- Sales number
- Serial number
- Expiry date
- Machine-dependent, maximum speed
- Maximum rotor cycles, maximum number of rotor cycles
- Current rotor cycles
- Remaining rotor cycles
- Minimum achievable temperature at maximum speed (only for centrifuges with cooling).

6.10.4.1.3 Error history

Personnel:

- Authorisation level 2
- Authorisation level 3

The *'Device information'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Error history'* menu is highlighted in blue.

2. ➤ Press *[Rotary, push button]*.

➤ Menu is displayed.

Warnings and errors that have occurred are displayed:

- Date of the warning or error
- Time of the warning or error
- Number of the warning or error

3. ➤ Rotate *[Rotary, push button]* until the desired error is highlighted in blue.

4. ➤ Press *[Rotary, push button]*.

➤ Error short text, error number and complete error text are displayed.

6.10.4.1.4 Operating hours

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'Device information'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *[Querying]* menu is highlighted in blue.

2. ➤ Press *[Rotary, push button]*.

➤ Menu is displayed.

The following information can be found in the menu:

- *'Current runtime'*
- *'Operating hours on the grid'*
- *'Total runtime'*

6.10.4.1.5 Data export to USB flash drive

'System log files' and 'Audit log files' can be exported to a USB flash drive from the menu.

The following file formats are supported for USB flash drives:

- FAT, FAT32, ext2, ext3, ext4

Personnel:

- Authorisation level 2
- Authorisation level 3

USB flash drive is inserted.

The 'Device information' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Data export to USB flash drive' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until the desired data for export is highlighted in blue.
4. ➤ Press [Rotary, push button].
 - Menu is displayed.
5. ➤ Rotate [Rotary, push button] until the desired time period for the export is highlighted in blue.
6. ➤ Press [Rotary, push button].
 - Menu with export requirements is displayed.
7. ➤ Press [Rotary, push button].
 - Export is started.

6.10.4.1.6 Service contact details

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'Device information' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Service contact details' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.

The following information can be found in the menu:

Service contact details:

- Phone:
- E-mail
- Note (information entered by service)

6.10.4.2 User settings

The user settings for authorisation levels 1 and 2 can be displayed only after a PIN has been entered by an authorisation level 3 user.

6.10.4.2.1 Starting point in Quick Adjust

A setting can be defined in the menu to determine whether the starting point in Quick Adjust is either the RPM/RCF or the time.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Starting point in Quick Adjust' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until the desired setting is selected.
4. ➤ Press the [BACK] button.
 - The settings are stored.

6.10.4.2.2 Runtime display

A setting can be defined in the menu to determine whether the runtime is displayed in 'enlarged' or 'classic' during the centrifugation run.

'enlarged':

The enlarged runtime display is shown during a centrifugation run or continuous centrifugation. In enlarged view, no changes can be made to the parameters during the centrifugation run.

'classic':

The classic runtime display is shown during a centrifugation run or during continuous centrifugation.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Runtime display' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until the desired setting is selected.
4. ➤ Press the [BACK] button.
 - The settings are stored.

6.10.4.2.3 Increment size for runtime settings

A setting can be defined in the menu to determine whether the increments when setting the runtime are 1 second or 15 seconds long.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The '*User settings*' submenu is selected in the settings menu.

1.  Rotate [*Rotary, push button*] until the '*Increment size for runtime setting*' menu is highlighted in blue.
2.  Press [*Rotary, push button*].
 - ➡ Menu is displayed.
3.  Rotate [*Rotary, push button*] until the desired setting is selected.
4.  Press the [*BACK*] button.
 - ➡ The settings are stored.

6.10.4.2.4 Opening the lid after the centrifugation run

The setting "Lid opens after centrifugation" is only available for Mikro 2.0.

A setting can be defined in the menu to specify whether the centrifuge lid opens automatically after the centrifugation run has ended.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The '*User settings*' submenu is selected in the settings menu.

1.  Rotate [*Rotary, push button*] until the '*Lid opens after centrifugation*' menu is highlighted in blue.
2.  Press [*Rotary, push button*].
 - ➡ Menu is displayed.
3.  Rotate [*Rotary, push button*] until the desired setting is selected.
4.  Press the [*BACK*] button.
 - ➡ The settings are stored.

6.10.4.2.5 Lid opens when switched on

A setting can be defined in the menu to specify whether the centrifuge lid opens automatically when the centrifuge is switched on.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The '*User settings*' submenu is selected in the settings menu.

1.  Rotate [*Rotary, push button*] until the '*Lid opens when switched on*' menu is highlighted in blue.
2.  Press [*Rotary, push button*].
 - ➡ Menu is displayed.

3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.6 Automatic start after lid closure

A setting can be defined in the menu to specify whether the centrifugation run starts automatically after the centrifuge lid has been closed.

The centrifugation run starts only when Quick Adjust or the program list is displayed.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Autom. start after lid closure'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.7 Service note

A setting can be defined in the menu to determine whether a service message should be displayed before the service interval expires. The service note appears after each centrifugation run.

The menu shows the number of days until the next service has to be carried out. If the service is overdue, a message appears.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Service note'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.8 Visual signal

A setting can be defined in the menu to determine whether a visual signal is displayed at the end of the centrifugation run. The visual signal is the gradual increase in screen brightness.

Opening the lid or pressing any button ends the signal.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Visual signal'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.9 Audible signal

A setting can be defined in the menu to determine whether a signal sounds after the centrifugation run is complete and in the event of warnings, and how loud the signals are.

'Audible signal at the end of centrifugation'

Three consecutive tones sound: "Low", "Medium" and "High".

- The tone sequence is repeated three times.
- This is followed by a pause of three seconds.
- The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

'Audible signal on warning'

Two consecutive tones sound: "Low" and "High".

- The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

'Audible signal on error'

The audible signal for errors cannot be configured. If an error occurs, the audible signal sounds at maximum volume.

- A high-pitched sound is emitted.
- The audible signal is repeated.

The signal is switched off by opening the lid or pressing any button.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Audible signal' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
 - 'Audible signal at the end of centrifugation'. The setting is highlighted in blue.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press *[Rotary, push button]*.
 - 'Audible signal on warning'. The setting is highlighted in blue.
5. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
6. ➤ Press *[Rotary, push button]*.
 - 'Audible signal volume'. The setting is highlighted in blue.
7. ➤ Rotate *[Rotary, push button]* until the desired audible signal volume is selected.
 - The audible signal is output during adjustment to allow the volume to be checked.

The volume setting governs the audible signal at the end of centrifugation and the audible signal for warnings.
8. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.10 Display animations

Display animations can be enabled and disabled from the menu.

Display animations are the transitions between different interfaces. If the setting is active, smooth transitions are displayed.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Display animations' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.11 Screen brightness

The screen brightness can be set in the menu.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Screen brightness' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired brightness is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.12 Energy saving mode

Energy saving mode can be set in the menu.

If the setting is active, the display switches off after the entered time if no button is pressed and no messages or signals appear.

Activate

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Energy saving mode' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until 'Yes' is selected.
4. ➤ Press *[Rotary, push button]*.
 - The time set for activation of energy-saving mode is displayed.
5. ➤ Rotate *[Rotary, push button]* until the desired time is set.

The time can be set in 30-minute increments up to 8 hours.
6. ➤ Press the *[BACK]* button.
 - The settings are stored.

Deactivate

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the 'Energy saving mode' menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until 'No' is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.13 Time

The time format can be selected and the time set in the menu.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Time' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until the desired time format setting is selected.
4. ➤ Press [Rotary, push button].
 - The next setting is highlighted in blue and displayed in bold.
5. ➤ Rotate [Rotary, push button] until the desired value for the hours is selected.
6. ➤ Press [Rotary, push button].
 - The next setting is highlighted in blue and displayed in bold.
7. ➤ Rotate [Rotary, push button] until the desired value for the minutes is selected.
8. ➤ Press the [BACK] button.
 - The settings are stored.

6.10.4.2.14 Date

The date format can be selected and the date set in the menu.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Date' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until the desired date format setting is selected.
4. ➤ Press [Rotary, push button].
 - The next setting is highlighted in blue and displayed in bold.
5. ➤ Rotate [Rotary, push button] until the desired value for the day or year is selected, depending on the date format set.
6. ➤ Press [Rotary, push button].
 - The next setting is highlighted in blue and displayed in bold.
7. ➤ Rotate [Rotary, push button] until the desired value for the month is selected.
8. ➤ Press [Rotary, push button].
 - The next setting is highlighted in blue and displayed in bold.

9. ➤ Rotate *[Rotary, push button]* until the desired value for the day or year is selected, depending on the date format set.
10. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.15 Temperature unit

The temperature unit setting is only available for the Mikro 2.0 R.

A setting can be defined in the menu to determine whether the temperature is displayed in degrees Celsius (°C) or degrees Fahrenheit (°F).

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Temperature unit'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired setting is selected.
4. ➤ Press the *[BACK]* button.
 - The settings are stored.

6.10.4.2.16 System language

The system language can be selected in the menu.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'System language'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - Menu is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired language is highlighted in blue.
4. ➤ Press *[Rotary, push button]*.
 - The language is selected.
5. ➤ Press the *[BACK]* button.
 - The settings are stored.

Changes take effect after a restart.

6.10.4.2.17 Authorisation level

Device protection requirements

The software of the MIKRO 2.0 and MIKRO 2.0 R includes appropriate measures to protect against unauthorised modification of parameters or system configurations due to (reasonably predictable) malicious third-party attempts.

If the device requires increased protection, the customer must provide additional measures to protect the device from unauthorised access from within the device's operating environment.

It is recommended that the MIKRO 2.0 and MIKRO 2.0 R be stored and used in restricted-access premises that limit device access to authenticated and authorised users.

Ensure that the device is intact before commissioning.

Some centrifuge functions cannot be performed by users with inadequate authorisation levels.

The authorisation level can be changed in the menu.

For authorisation level 1

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Authorisation level' menu is highlighted in blue.
2. ➤ Press [Rotary, push button].
 - Menu is displayed.
3. ➤ Rotate [Rotary, push button] until Authorisation Level 1 is highlighted in blue.
4. ➤ Press [CONTEXT BUTTON, TOGGLE BUTTON].
 - Context menu is displayed.
5. ➤ Rotate [Rotary, push button] until the desired program, which should be retrievable at authorisation level 1, is highlighted in blue.
6. ➤ Press [Rotary, push button].
 - The program is selected.
 - Any number of programs can be selected.
7. ➤ Press the [BACK] button.
 - The permitted programs are saved.
8. ➤ Press the [BACK] button.
 - The settings are stored.

For authorisation levels 2 and 3

Personnel:

- Authorisation level 2
- Authorisation level 3

The 'User settings' submenu is selected in the settings menu.

1. ➤ Rotate [Rotary, push button] until the 'Authorisation level' menu is highlighted in blue.

2.  Press *[Rotary, push button]*.
 - ➡ Menu is displayed.
3.  Rotate *[Rotary, push button]* until the authorisation level is highlighted in blue.
4.  Press *[Rotary, push button]*.
 - ➡ The authorisation level is selected.
5.  Press the *[BACK]* button.
 - ➡ The settings are stored.

6.10.4.2.18 Changing the PIN

The PIN for users with authorisation level 3 can be changed in the menu.

Personnel:

- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1.  Rotate *[Rotary, push button]* until the *'Changing the PIN'* menu is highlighted in blue.
2.  Press *[Rotary, push button]*.
 - ➡ The menu for entering the old PIN is displayed.
3.  Rotate *[Rotary, push button]* to select the desired number, then press *[Rotary, push button]* to enter the number.

To delete a number, rotate [Rotary, push button] until the  symbol is highlighted in blue, then press [Rotary, push button].

Repeat the process until all 8 numbers have been entered.

- ➡ ✓ symbol is highlighted in blue.

4.  Press *[Rotary, push button]*.
 - ➡ The menu for entering the new PIN is displayed.
5.  Rotate *[Rotary, push button]* to select the desired number, then press *[Rotary, push button]* to enter the number.

To delete a number, rotate [Rotary, push button] until the  symbol is highlighted in blue, then press [Rotary, push button].

Repeat the process until all 8 numbers have been entered.

- ➡ ✓ symbol is highlighted in blue.

6.  Press *[Rotary, push button]*.
 - ➡ The menu for confirming the new PIN is displayed.

7. ➤ Rotate *[Rotary, push button]* to select the desired number, then press *[Rotary, push button]* to enter the number.

*To delete a number, rotate *[Rotary, push button]* until the  symbol is highlighted in blue, then press *[Rotary, push button]*.*

Repeat the process until all 8 numbers have been entered.

- ✓ symbol is highlighted in blue.

8. ➤ Press *[Rotary, push button]*.

- The settings are stored.

Information line indicates successful execution of the action.

If you forget your PIN, you can reset it using the advanced settings and an authentication file available from Service.

6.10.4.2.19 Overwriting rotor info

In the menu, information from a previously used rotor can be replaced by information from the newly inserted rotor.

The function is available when memory locations for the rotor information are full. This is signalled by warning message 2.2.300.018.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The *'User settings'* submenu is selected in the settings menu.

1. ➤ Rotate *[Rotary, push button]* until the *'Overwriting rotor info'* menu is highlighted in blue.
2. ➤ Press *[Rotary, push button]*.
 - The list of rotors is displayed.
3. ➤ Rotate *[Rotary, push button]* until the desired rotor is highlighted in blue.
4. ➤ Press *[Rotary, push button]*.
 - Confirmation query appears.
5. ➤ Rotate *[Rotary, push button]* until symbol  is highlighted in blue.
6. ➤ Press *[Rotary, push button]*.
 - The rotor information is overwritten.

Information line indicates successful execution of the action.
7. ➤ Press the *[BACK]* button.
 - The submenu is exited.

6.10.4.2.20 Default settings

The menu allows all settings to be reset to their default settings.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The '*User settings*' submenu is selected in the settings menu.

1. ➤ Rotate [*Rotary, push button*] until the '*Default settings*' menu is highlighted in blue.
2. ➤ Press [*Rotary, push button*].
 - Menu is displayed.
 - '*Continue*' is highlighted in blue.
3. ➤ Press [*Rotary, push button*].
 - Confirmation query appears.
4. ➤ Rotate [*Rotary, push button*] until symbol ☑ is highlighted in blue.
5. ➤ Press [*Rotary, push button*].
 - User settings are reset to their defaults.
 - Information line indicates successful execution of the action.
6. ➤ Press the [*BACK*] button.
 - The submenu is exited.

6.10.4.3 Service settings and Advanced settings

The '*Service settings*' and '*Advanced settings*' submenus cannot be accessed by users with authorisation levels 1, 2 or 3.

These submenus are reserved for customer service and so are not described in the operating instructions.

To exit the authentication query, press [*Rotary, push button*] when the ☒ symbol is highlighted in blue.

If the ☒ symbol is not highlighted in blue, first rotate [*Rotary, push button*] so that the ☒ symbol is highlighted in blue.

6.10.4.4 Open source licences

The open source licences used can be displayed in the menu.

Personnel:

- Authorisation level 1
- Authorisation level 2
- Authorisation level 3

The '*Open source licences*' submenu is selected in the settings menu.

1. ➤ Rotate [*Rotary, push button*] to scroll up and down through the open source licences.
2. ➤ Press the [*BACK*] button.
 - The submenu is exited.

7 Maintenance, cleaning and care

! WARNING

Risk of electric shock due to maintenance and servicing work on live device.

- Disconnect the device from the mains before carrying out repairs and maintenance.

DANGER

Risk of contamination

Inadequate cleaning or failure to observe the cleaning instructions can lead to contamination risks.

- Observe national and local regulations on safety and accident prevention.
- Observe cleaning instructions.
- Wear personal protective equipment when cleaning the device.
- Comply with laboratory regulations (e.g. TRBAs, the German Protection against Infection Act, hygiene plan) for handling biological agents.

Instructions for cleaning and disinfection

- The device and its accessories must not be cleaned in dishwashers.
- Only perform hand cleaning and liquid disinfection.
- The water temperature must not exceed 25 °C.
- To prevent any corrosion due to use of detergents or disinfectants, it is essential to follow the special application instructions provided by the manufacturers of the detergent or disinfectant.

Disinfectant

- Surface disinfectant (not disinfectant for hands or instruments)
- Ethanol as the sole active substance. Do not use an ethanol-propanol mixture to disinfect the viewing window in the lid of the device.
- Concentration is not less than 30%
- pH: 6 – 8
- Non-corrosive

7.1 Cleaning

Maintenance work	Interval in operating hours (h)	To be performed by
Cleaning the device	8	Trained person
Cleaning biosafety systems	100	Trained person
Cleaning accessories	100	Trained person

Cleaning the device

1. Open the lid.
2. Switch off the device and disconnect it from the power supply.
3. Remove accessories.
4. Clean the centrifuge housing and the centrifuging chamber with soap or a mild detergent and a damp cloth.

5. ➤ Remove any detergent residues with a damp cloth after using detergents.
6. ➤ The surfaces must be dried immediately after cleaning.
7. ➤ Dry the centrifuging chamber with an absorbent cloth if condensation forms.

Cleaning biosafety systems

1. ➤ Clean the biosafety system using the detergent and a damp cloth.
2. ➤ Remove any detergent residues with a damp cloth after using detergents.
3. ➤ Dry the accessories immediately after cleaning using a lint-free cloth and oil-free compressed air. Dry all cavities completely using oil-free compressed air.

Cleaning accessories

1. ➤ Clean the accessories using the detergent and a damp cloth.
2. ➤ Remove any detergent residues with a damp cloth after using detergents.
3. ➤ Dry the accessories immediately after cleaning using a lint-free cloth and oil-free compressed air. Dry all cavities completely using oil-free compressed air.

7.2 Disinfection

Disinfection must always be preceded by cleaning of the components concerned.

Disinfectant concentration and application time according to the manufacturer's instructions.

Maintenance work	Interval in operating hours (h)	To be performed by
Disinfecting the device	if required	Trained person
Disinfecting the accessories	if required	Trained person
Autoclaving	if required	Trained person

Disinfecting the device

CAUTION

Risk of electric shock

- Liquids that penetrate the device may cause electric shocks.
- The device must be protected from external contact with liquids.
 - Do not pour any liquids inside the device.
 - The original transport packaging must be used when transporting the device.

1. ➤ Open the lid.
2. ➤ Switch off the device and disconnect it from the power supply.
3. ➤ Remove accessories.
4. ➤ Clean the housing and centrifuging chamber using disinfectant.
5. ➤ Remove any disinfectant residues with a damp cloth after using disinfectants.
6. ➤ The surfaces must be dried immediately after cleaning.

Disinfecting the accessories

1. ➤ Disinfect the accessories using the disinfectant.
2. ➤ Wet all cavities with bubble-free disinfectant.
3. ➤ Remove the disinfectant residues or leave them to dry after using disinfectants.

Autoclaving

The following accessories can be autoclaved at 134 °C / 273 °F, 2 bar (10 min) up to 10 times:

- Aluminium angle rotors
- Lid

Before autoclaving, the rotor lid must be removed and the seals removed from the rotor.

No statement can be made about the resulting degree of sterility.

Autoclaving accelerates the ageing of materials. It may cause changes to colours.

1. ➤ Remove the lid of the rotors before autoclaving.
2. ➤ After autoclaving, visually check rotors and accessories for damage.
3. ➤ Replace damaged parts.
4. ➤ Replace the sealing ring if it is cracked, brittle or worn.
5. ➤ For biosafety systems, replace the sealing rings after autoclaving.

7.3 Maintenance

WARNING

There is a risk of damage to the cooling circuit.
The cooling circuit must not be damaged.

Maintenance work	Interval in operating hours (h)	To be performed by
Checking the accessories	100	Qualified mechanics
Checking the biosafety system	100	Qualified mechanics
Centrifuging chamber damage inspection	Annually	Qualified mechanics
Accessories with a limited service life	if required	Qualified mechanics
Replacing centrifuge tubes	if required	Qualified mechanics

Checking the accessories

1. ➤ The accessories are to be checked for wear and corrosion damage.
 - Remove accessories from operation if they show signs of wear.

2. ➤ Check that the rotor is firmly seated.

Checking the biosafety system

1. ➤ Visually check all parts of the biosafety system for damage.
2. ➤ Check the correct installation position of the sealing ring(s) of the biosafety system.
3. ➤ Replace the damaged parts of the biosafety system.
4. ➤ Replace the sealing ring if it is cracked, brittle or worn. Replace lids with fixed sealing rings as a complete unit.

Centrifuging chamber damage inspection

- Check the centrifuging chamber for damage.

Accessories with a limited service life

Do not use accessories if the maximum number of running cycles or the expiry date has been reached.

- The maximum permissible number of running cycles or the expiry date must be complied with.
- The centrifuge is equipped with a cycle counter.

Replacing centrifuge tubes

CAUTION

Risk of injury

Glass breakage and contaminated liquids in the centrifuge.

- Wear personal protective equipment.
- Stop the centrifuge immediately and disconnect it from the power supply.
- No contact with broken glass or spilled substances.
- Clean the area in accordance with the safety and disposal regulations, watching out for sharp objects and biological material.
- Observe national and local environmental protection and disposal regulations to ensure proper disposal.

Broken parts of the tube, glass fragments and spilled centrifuge material must be removed completely in the event of leakage or if a centrifuge tube breaks. Glass splinters that are not removed will cause further glass breakage.

The rubber inserts and the plastic sleeves of the rotors must be replaced after a glass breakage.

Disinfection must be carried out if the material is infectious.

Rotor sealing ring replacement

Remove inner seal and reinstall

Personnel:

- Trained user

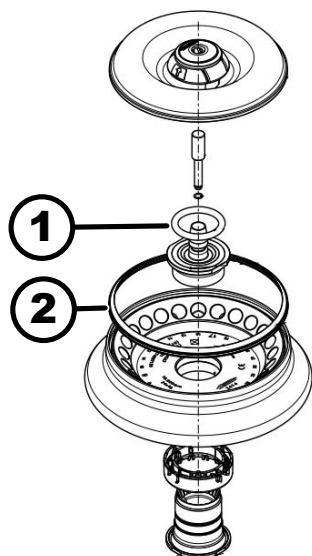
Fig. 24: Rotor sealing ring positions

- 1 Inner seal
- 2 Outer seal: Lip seal

1. ➤ Remove the inner seal from the groove.
Do not use pointed or sharp objects.
2. ➤ Check the groove for cleanliness and damage.
3. ➤ Check the seal for damage before inserting it.
4. ➤ Insert the seal into the groove.

Remove outer seal (lip seal) and reinstall

Personnel:

- Trained user

1. ➤ Pull the lip seal out of the groove.
2. ➤ Check the groove for cleanliness and damage.
3. ➤ Check the lip seal for damage before inserting it.
4. ➤ Align the lips of the lip seal upwards and insert smoothly and without creases into the groove.

8 Troubleshooting

8.1 Fault description

Personnel:

- Trained person

In the event of faults that are not listed, contact customer service. Specify the centrifuge type and serial number from the centrifuge rating plate.

Error code number	Cause	Remedy
1.1.001.000 - 1.1.002.004	Error in the tachometer system	Wait until the centrifuge comes to a standstill. Then open the lid. Then switch off the centrifuge. Rotate the rotor by hand and switch on the centrifuge. If the error message cannot be deleted, contact Service.

Error code number	Cause	Remedy
1.1.004.001 - 1.1.004.005	Lid closure error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.005.000 - 1.1.005.003	Error due to overspeed	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.006.000	Lid closure time-out	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.007.001, 1.1.007.002	Lid closure control error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.010.001 - 1.1.010.005	Error in rotor detection	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.011.001	Power supply interruption	Clear the error by performing a power reset.
1.1.011.002	Mains voltage interruption detected at standstill	Clear the error by performing a power reset. If the fault occurs again, undertake mains voltage stabilising measures. If the mains voltage is stable, contact Service.
1.1.012.001, 1.1.012.002	Invalid machine version	Contact service.
1.1.012.003	Invalid machine component DFC	Contact service.
1.1.012.004	Invalid machine component CFC	Contact service.
1.1.012.005	Invalid machine component DFC	Contact service.
1.1.012.006	Invalid machine component CFC	Contact service.
1.1.014.002	Target speed exceeded	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.014.003	Actual speed too low	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.025.001	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.025.002 - 1.1.025.006	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.026.000 - 1.1.026.050	Sensor defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.

Error code number	Cause	Remedy
1.1.027.000, 1.1.027.001	System test Centrifuge control unit error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.028.001 - 1.1.028.007	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.030.001, 1.1.030.002	Centrifuge control unit error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.031.000 - 1.1.049.003	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.051.000 - 1.1.055.000	Error in the cooling system	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.057.000	Temperature control failed	The manufacturer recommends compliance with the temperature control specifications. If this occurs repeatedly, contact Service.
1.1.058.000 - 1.1.058.007	Error in the cooling system	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.060.000	Drive frequency converter error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.060.001	DFC error due to closure	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.061.000, 1.1.061.001	DFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.062.000, 1.1.062.001	CFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.063.000 - 1.1.063.003	DFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.063.004	CFC error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.070.000, 1.1.070.001	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.

Error code number	Cause	Remedy
1.1.071.000, 1.1.071.001	[STOP, OPEN] button defective	The [STOP, OPEN] button was identified as defective after switching on the centrifuges. Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.086.000 - 1.1.086.012	Internal error in the centrifuge control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.091.000	Imbalance sensor X-axis defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.091.001, 1.1.091.002	Imbalance sensor X-axis offset error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.092.000	Imbalance sensor Y-axis defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.092.001, 1.1.092.002	Imbalance sensor Y-axis offset error	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.093.000	Temperature sensor in imbalance sensor defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.096.000 - 1.1.096.004	Error in the tachometer system	Delete the error message: Wait until the centrifuge comes to a standstill. Then open the lid. Then switch off the centrifuge. Rotate the rotor by hand and switch on the centrifuge. If the error message cannot be deleted, contact Service.
1.1.097.000, 1.1.098.000	Sensor defective	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.521.001	Communication error, control unit and drive	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
1.1.521.003	NAK for communication between control unit and drive	Contact service.
1.1.521.004	No communication between control unit and drive	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.

Error code number	Cause	Remedy
1.1.521.005, 1.1.521.006	Communication error drive control unit	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.400.001	[STOP, OPEN] button malfunction detected	Contact service.
2.1.420.001	Centrifuge control unit type error	Switch off the device. Contact service.
2.1.501.001	Centrifuge control unit illogical information	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.541.001	Power supply interruption	Clear the error by performing a power reset.
2.1.551.001	Error message logging	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.600.001	Error in the application	This error message can only be deleted by performing a power reset.
2.1.620.000	ServiceAuth file incorrect	Clear the error by performing a power reset.
2.1.800.002 - 2.1.800.010	Error message logging	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.002	Drive is not ready	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.003	Function test not possible	Clear the error by performing a power reset. If this occurs repeatedly, contact Service.
2.1.901.004	Centrifuge control unit version error	Contact service.
2.1.901.005	Centrifuge control unit configuration error	Contact service.

8.2 Warning descriptions

Personnel:

- Trained person

In the event of faults that are not listed, contact customer service. Specify the centrifuge type and serial number from the centrifuge rating plate.

Warning code number	Cause	Remedy
1.2.003.001	The centrifuge has detected an imbalance.	Please open the lid when the centrifuge is at a standstill. Check loading. Ensure even weight distribution.
1.2.011.000	A power failure occurred during centrifugation.	Open the lid when the centrifuge is at a standstill. The manufacturer recommends checking the result of the centrifugation.
2.2.300.001	The centrifuge has detected a rotor change.	Acknowledgement of the warning only possible at standstill. The centrifuge will check or adjust the current centrifugation parameters for the current rotor.
2.2.300.002	Starting of centrifugation without a rotor.	Acknowledgement of the warning only possible at standstill. The centrifuge has detected that centrifugation has been started without a rotor. Only start centrifugation with a rotor.
2.2.300.003	The permissible number of cycles for the rotor currently in use has been reached.	The rotor currently in use must be replaced immediately with a new rotor. We warn against further use. Further use only at your own risk.
2.2.300.004	The permissible number of cycles for the rotor currently in use has been reached.	Use of the current rotor poses a safety risk. Use only at your own risk.
2.2.300.005	The centrifuge has detected a rotor change.	The centrifuge will check or adjust the current centrifugation parameters for the current rotor.
2.2.300.010, 2.2.300.011	The permitted number of centrifugations has been reached in spite of a faulty rotor identification system.	Further centrifugation is not possible. Contact service.
2.2.300.012	The rotor identification system (tag) is faulty or no rotor is fitted.	The number of centrifugation cycles is limited in this state. It is strongly recommended that you contact Service if the rotor identification system malfunctions. If no rotor is fitted, the start of centrifugation is terminated with a warning message.
2.2.300.013	The permissible operating period for the rotor has been exceeded.	The rotor currently in use must be replaced immediately with a new rotor. We warn against further use. Further use only at your own risk.

Warning code number	Cause	Remedy
2.2.300.014	The rotor identification system (reader) is faulty.	The number of centrifugation cycles is limited in this state. Contact service.
2.2.300.015	The rotor identification system reports an error.	The number of centrifugation cycles is limited in this state. Contact service. More information on the number of centrifugations at the start of this.
2.2.300.016	The current rotor is not approved for this centrifuge. It is not possible to start centrifugation.	Approved rotors can be found in the operating instructions. Contact service.
2.2.300.017	The current rotor does not comply with the Hettich standard.	It is not possible to start centrifugation. Please remove this rotor from the centrifuge immediately. Contact service.
2.2.300.018	The memory locations for rotor information are full. It is currently not possible to start centrifugation.	Select an entry in "Overwrite rotor info" that may be overwritten with the data of the current rotor. This action is only possible when the new rotor is in the centrifuge.
2.2.300.019, 2.2.300.020	The permitted number of centrifugations has been reached in spite of a faulty rotor identification system.	The manufacturer recommends reactivating the centrifugation allocation. Contact service.
2.2.310.001, 2.2.310.002	The user has cancelled the current centrifugation by pressing the <i>[STOP, OPEN]</i> button.	Acknowledgement of the warning only possible at standstill. The manufacturer recommends checking the result of the centrifugation.
2.2.320.001	User has cancelled the data export.	Remove USB flash drive or restart data export.
2.2.340.001	The desired temperature is above the current temperature. This centrifuge can only cool.	To achieve a higher temperature than the current one, carry out a standard centrifugation.
2.2.550.013	The update package does not contain a more up-to-date software version for the centrifuge.	The software update is cancelled.
2.2.550.014	The software version of the HBI update file was not accepted.	Check the version of the software update used.
2.2.550.015	The ZS software version has not been updated.	Repeat software update.

Warning code number	Cause	Remedy
2.2.560.001	Failed to mount the USB flash drive.	Use another USB flash drive. If this occurs repeatedly, contact Service.
2.2.560.002	Unmounting the USB flash drive failed.	Remove the USB flash drive and restart the centrifuge. If this occurs repeatedly, contact Service.
2.2.560.003	The system reports: "No USB flash drive available".	Remove the USB flash drive and insert it again. Use a different USB flash drive if the warning message appears again. If there is a permanent warning message, contact Service.
2.2.560.004	Insufficient storage space on USB flash drive for data export.	Use a USB flash drive with sufficient free storage space. Recommendation: at least 8 GB
2.2.560.005	Failed to write to USB flash drive. Data may not have been completely written to the USB flash drive.	Check USB flash drive and restart export.
2.2.560.006	The file system of the USB flash drive does not correspond to the system specifications. Data storage is only possible with the file systems: FAT, FAT32, ext2, ext3, ext4.	Use a USB flash drive with a supported file system.
2.2.560.007	Deletion of a file on the USB flash drive has failed. Copying process was cancelled.	Clean up the USB flash drive and export again.
2.2.560.008	Remounting the USB flash drive has failed. Files may not have been copied completely.	Check USB flash drive and perform the export again.
2.2.570.001	Next service imminent.	Contact service.
2.2.600.002	A defective database was detected and restored.	Acknowledge warning message. If this occurs repeatedly, contact Service.
2.2.600.003	The database restore has failed. The database has been recreated. Programs and settings are no longer available.	Contact service.
2.2.600.004	Error when deleting the system logs.	Perform "Reset to delivery status" again. If this occurs repeatedly, contact Service.

Warning code number	Cause	Remedy
2.2.600.010	A problem with the date or time was detected when the centrifuge was started.	Please correct the date or time in centrifuge. If this occurs repeatedly, contact Service.
2.2.600.011	The battery of the internal clock is discharged. The service request was activated out of schedule.	To replace the internal battery, contact Service.
2.2.600.021	Machine version rejected.	Limited functionality, no centrifugation possible. Contact service.
2.2.600.040	Balancing of the motor and DFC is mandatory.	Centrifugation is only possible after balancing. Contact service.

8.3 Performing a MAINS RESET

1. ➤ Set the mains switch to [0].
2. ➤ Wait 10 seconds.
3. ➤ Set the mains switch to [I].

8.4 Emergency release

The lid cannot be unlocked by the motor in the event of a power failure. Emergency unlocking by hand must be performed.

! WARNING

Risk of electric shock due to maintenance and servicing work on live device.

- Disconnect the device from the mains before carrying out repairs and maintenance.


WARNING

Risk of cuts and crushing

Risk of injury due to moving rotor.

- Do not open the lid until the rotor has come to a complete stop.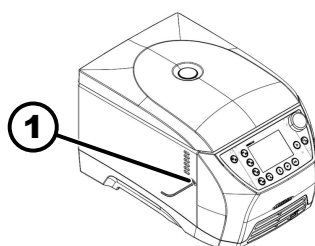

1 Hole

Personnel:

- Trained user

1. ➤ Look through the window in the lid to ensure that the rotor is stationary.
2. ➤ Insert an Allen key horizontally into the hole (1).
3. ➤ Rotate the Allen key anticlockwise until the lid opens.

8.5 Replacing the mains input fuse

! WARNING

Risk of electric shock due to maintenance and servicing work on live device.


- Disconnect the device from the mains before carrying out repairs and maintenance.

Personnel:

- Trained user

The mains fuses are located next to the mains switch.

The mains switch is in switch position [0]

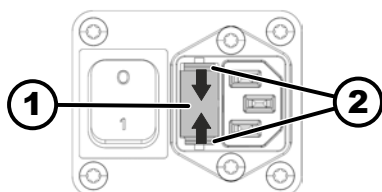- 1 Fuse holder
- 2 Snap lock

1. ➤ Disconnect the mains cable from the device plug.
2. ➤ Press the snap locks (2) against the fuse holder (1) and pull them out.

3. ➤ Replace the defective mains input fuses.

Only use fuses with the nominal value specified for the type: see the table below.

4. ➤ Push in the fuse holder (1) until the snap lock engages.

5. ➤ Reconnect the device to the mains.

Model	Type	Fuse	Order no.
MIKRO 2.0	2404	T8 AH/250V	E738
	2407		
MIKRO 2.0 R	2406	T8 AH/250V	E738
	2408		

9 Spare parts

Only use original spare parts from the manufacturer, approved accessories and wear parts.

10 Disposal

10.1 General instructions

The device can be disposed of via the manufacturer.

A Return Material Authorisation (RMA) form must always be requested for a return.

If necessary, contact the Technical Service Department of the manufacturer:

- **Andreas Hettich GmbH**
- Föhrenstraße 12
- 78532 Tuttlingen, Germany
- Phone: +49 7461 705 1400
- E-Mail: service@hettichlab.com

The cooling circuit must only be dismantled by qualified personnel.

No sensitive data (i.e. data requiring protection, such as patient information) is stored on the MIKRO 2.0. A factory reset is recommended before decommissioning the device in order to safely remove program data/configurations from the system. ➔ Chapter 6.10.4.2.20 'Default settings' on page 67

! WARNING

Fire and explosion hazard due to flammable refrigerant (R290, safety class A3 according to ISO 817 and ASHRAE 34).

The refrigerant must be evacuated before dismantling the refrigerant circuit.

- Keep refrigerant away from ignition sources.

! WARNING

Risk of soiling and contamination

Damage to the environment and health due to improper disposal.

- Observe national and local environmental protection and disposal regulations to ensure proper disposal or recycling. Separate metals, non-metals, composite materials and auxiliary materials by type and dispose of them in an eco-friendly manner.
- The device may only be disassembled and disposed of by trained and authorised service technicians.

The device is intended for the commercial sector ("Business to Business" - B2B).

According to Directive 2012/19/EU, the devices may no longer be disposed of with household waste.

The devices are assigned to the following groups according to the Stiftung Elektro-Altgeräte Register (EAR (German foundation under civil law)):

■ Group 5 (small devices)

The crossed-out wheellie bin symbol indicates that the device must not be disposed of with household waste. Regulations governing disposal of such devices may differ in individual countries. If necessary, contact the supplier.

Fig. 25: Household waste ban

11 Free and Open Source Software

This product contains free and open source software. See the enclosed DVD "Licenses and Sourcecode" for the the applicable license texts, the source code, copyright notices for the individual components, any disclaimers, permission notes and other information about the individual components. The DVD contains the following folders:

File	Content
File u-boot-*.*	Folders with the license text of the uboot bootloader components
gelin2-19.10-imx6-x86_64.tar.*	Complete Linux packet (GELin BSP) incl. license info and source code
HettCube-GELin_KernelV5.4.53_Lizenzinformationen.txt	Links to the license info of the Linux Kernel V5.4.53
hettcube-gelin2-19.10_BoM.csv	List of all GELin components with detailed information to the licenses
hettcube-gelin2-19.10_license-information.*	License text files (txt and HTML)
Release.Kernel.linux-5.4.53-ge-20.08.for.exceet_imx6dl_ts_hmi-hettich.zip	Linux Kernel V5.4.33 (adapted for the use with the HMI electronic board)

Please note: Use of free and open source software components listed on the DVD files mentioned in the table above are governed by the terms of the licenses of the respective copyright holder as described in the folders on the DVD. Some of the licenses contain liability and warranty disclaimers and are provided "as is" without warranty of any kind. For more details and additional conditions for using the software, please read the attached licenses.

License Terms and Conditions regarding LGPL components:

For those parts of the HETTICH software that are licenses under

- the "GNU Library General Public License" or the "GNU Lesser General Public License" (LGPL) version 2.1 and earlier ("work") the customer is expressly allowed to make modifications of the work for the customer's own use and to reverse engineer for debugging such modifications;
- the "GNU Lesser General Public License" (LGPL) version 3 and later ("work") it is expressly allowed to make modifications of the portions of the Library contained in the work and reverse engineer for debugging such modifications.

Further rights of the customer under applicable copyright law are not restricted. Please note, that those parts of the software which are under free and open source licenses are distributed free of charge.

Installation Information for modified components:

Modified software components, esp. those under the GNU Lesser General Public License version (LGPL), can be installed by copying the new software component to the storage device contained within the product.

**CAUTION**

CHANGING THE SOFTWARE OF THE PRODUCT MAY RESULT IN INJURY TO LIFE, LIMB, OR HEALTH! ALSO IT CAN IMPACT THE CERTIFICATION OF THE PRODUCT AND RESULT IN A NON-LEGAL USE OF THE DEVICE. CHANGES CAN ALSO AFFECT THE WARRANTY.

Special License Information

The enclosed DVD "HettCube Licenses and Sourcecode" is part of the present documentation. We would like to point out the following separately:

OpenSSL

This product includes software developed by the OpenSSL Project for use in the OpenSSL Toolkit (<http://www.openssl.org>).

Eric Young

This product includes cryptographic software written by Eric Young (ey@cryptsoft.com).

Fontconfig

This product includes software under the fontconfig license.

- Copyright © 2000,2001,2002,2003,2004,2006,2007 Keith Packard
- Copyright © 2005 Patrick Lam
- Copyright © 2009 Roozbeh Pournader
- Copyright © 2008,2009 Red Hat, Inc.
- Copyright © 2008 Danilo Šegan
- Copyright © 2012 Google, Inc.

Permission to use, copy, modify, distribute, and sell this software and its documentation for any purpose is hereby granted without fee, provided that the above copyright notice appear in all copies and that both that copyright notice and this permission notice appear in supporting documentation, and that the name of the author(s) not be used in advertising or publicity pertaining to distribution of the software without specific, written prior permission. The authors make no representations about the suitability of this software for any purpose. It is provided "as is" without express or implied warranty.

THE AUTHOR(S) DISCLAIMS ALL WARRANTIES WITH REGARD TO THIS SOFTWARE, INCLUDING ALL IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS, IN NO EVENT SHALL THE AUTHOR(S) BE LIABLE FOR ANY SPECIAL, INDIRECT OR CONSEQUENTIAL DAMAGES OR ANY DAMAGES WHATSOEVER RESULTING FROM LOSS OF USE, DATA OR PROFITS, WHETHER IN AN ACTION OF CONTRACT, NEGLIGENCE OR OTHER TORTIOUS ACTION, ARISING OUT OF OR IN CONNECTION WITH THE USE OR PERFORMANCE OF THIS SOFTWARE.

The FreeType Project License

This product includes software based in part of the work of the FreeType Team (<https://freetype.org/>).

For further information (including license texts and copyright notices) see the enclosed DVD.

12 Glossary

Autoclaving	Autoclaving is a process for sterilising materials and objects using pressurised saturated steam in an autoclave. An autoclave is a closed device capable of generating high pressure and high temperatures to kill microorganisms such as bacteria, viruses, fungi and spores.
Brake level	Determines how quickly the centrifuge is decelerated from maximum speed to a complete standstill, influenced by various levels from gentle to fast.
Centrifugation	The process of applying centrifugal force to separate components of a solution based on their density.
Centrifugation allocation	Allocation (number) of approved centrifugations that the customer may perform despite a malfunction in the rotor identification system.
Commissioning	The process of setting up and starting the centrifuge before first use or after maintenance.
General information about the glossary	This glossary provides a comprehensive overview of key terms that are important when using the centrifuge and helps laboratory personnel to better understand the documentation and operating instructions for the centrifuge.
Lid	The part of the centrifuge that remains closed during operation to ensure safety and protect the integrity of the samples.
PIN	Personal identification number used to access secure centrifuge functions.
Power reset	A procedure for resetting the electrical supply to the centrifuge in order to correct system errors or restart the system.
Ramp-up level	Defines the rate of acceleration of the centrifuge up to the target speed. Different levels enable gentle or rapid acceleration depending on sample sensitivity.
RCF (Relative centrifugal force)	relative centrifugal force The force acting on the sample during centrifugation, measured as a multiple of the acceleration due to gravity (g).
Residual current circuit breaker (RCD)	The residual current circuit breaker interrupts the power supply in the event of a residual current and thus protects against electric shocks and fires.
Risk of contamination	Risk of contamination of the samples or the working environment, in particular by pathogenic organisms or chemical substances.
Rotor	The rotating part of the centrifuge in which the sample containers are inserted. Rotors can vary depending on the application.
Safety area	The area around the centrifuge that must be free of obstacles to ensure safety during operation.
UDI (Unique Device Identification)	A worldwide system for the unique identification of medical devices over their entire service life.

13 Index

A

Accessories

autoclaving.	71
biosafety systems.	70
checking.	71
disinfection.	71

Authorisation levels

Authorisations.	8
----------------------	---

Autoclaving.	71
-------------------	----

B

Biosafety system

checking.	72
----------------	----

Brake level.	30
-------------------	----

C

Centrifugation

Short-term centrifugation.	39
with higher substance density.	29
with time preselection.	38

Centrifuge tubes

filling.	27
replacing.	72

Certifications.	14
----------------------	----

Cleaning

biosafety systems.	70
-------------------------	----

Connecting the centrifuge.	22
---------------------------------	----

Continuous operation.	37
----------------------------	----

D

Device

biosafety systems.	69
disinfection.	70

Device information

Data export to USB flash drive.	56
Error history.	55
Querying.	54, 55
Rotor information.	54
Service contact details.	56

Disposal.	82
----------------	----

E

Error messages.	73
----------------------	----

F

Filling.	27
---------------	----

G

General safety instructions.	10
-----------------------------------	----

I

Information lines.	33
-------------------------	----

L

Labels

on the device.	16
on the packaging.	15

Lid

the lid.	25
---------------	----

Loading.	27
Logos.	14

M

MAINS RESET.	81
-------------------	----

Menus

Main menu.	35
Program menu.	37
Quick Adjust menu.	36
Settings menu.	37

O

Opening

a program link.	44
changing.	45
copying.	50
deletion.	52
Pre-cooling.	43
renaming.	51

Operator obligations

Obligations of the operator.	9
-----------------------------------	---

P

Program

retrieving.	45
------------------	----

Program favourite

Creating favourites.	47
deletion.	49

Program favourites

retrieving.	48
------------------	----

Q

Quick Adjust menu.	41
-------------------------	----

R

Ramp-up level.	30
---------------------	----

Rating plate.	13
--------------------	----

Relative centrifugal force

RCF.	29
-----------	----

Removing

Symbols.	28
the rotor.	28

Return.	16
--------------	----

Rotor

inserting.	25
removing.	25

Rotor lid

fitting.	26
removal.	26

Runtime.	30
---------------	----

Runtime starting time

changing.	42, 47
----------------	--------

S

Safety instructions.	10
---------------------------	----

Scope of delivery.	18
-------------------------	----

Setting up the centrifuge.	20
---------------------------------	----

Settings menu

Advanced settings.	68
Device information.	54

Open source licences.	68
Service settings.	68
User settings.	57
Speed.	42, 46
Storage conditions.	19
Switching off.	24
Switching the unit on.	24
Switching on for the first time.	22
Symbols.	6, 32
T	
Transport condition.	18
Transport lock	
fastening.	20
removal.	20
Troubleshooting.	73
U	
Unpacking.	20
User interface	
Keypad - Delete input.	35
Keypad - letters, numbers.	34, 35
Keypad - space bar.	34
Keypad - special characters.	35
Keypad - Upper case letters.	34
User settings	
Audible signal.	60
Authorisation level.	65
Automatic start after closing the lid.	59
Changing the PIN.	66
Date.	63
Default settings.	67
Display animations.	61
Energy saving mode.	62
Increment size for runtime settings.	58
Lid opens after centrifugation.	58
Lid opens when switched on.	58
Overwriting rotor info.	67
Runtime display.	57
Screen brightness.	61
Service note.	59
Starting point in Quick Adjust.	57
System language.	64
Temperature unit.	64
Time.	62
Visual signal.	60

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File	Content
File u-boot-*.*	Folders with the license text of the uboot bootloader components
gelin2-19.10-imx6-x86_64.tar.*	Complete Linux packet (GELin BSP) incl. license info and source code
HettCube-GELin_KernelV5.4.53_Lizenzinformationen.txt	Links to the license info of the Linux Kernel V5.4.53
hettcube-gelin2-19.10_BoM.csv	List of all GELin components with detailed information to the licenses

File	Content
hettcube-gelin2-19.10_license-information.*	License text files (txt and HTML)
Release.Kernel.linux-5.4.53-ge-20.08.for.exceet_imx6dl_ts_hmi-hettich.zip	Linux Kernel V5.4.33 (adapted for the use with the HMI electronic board)

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Eric Young

This product includes cryptographic software written by Eric Young (eay@cryptsoft.com).

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11 Free and Open Source Software

This product contains free and open source software. See the enclosed DVD "Licenses and Sourcecode" for the applicable license texts, the source code, copyright notices for the individual components, any disclaimers, permission notes and other information about the individual components. The DVD contains the following folders:

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Eric Young

This product includes cryptographic software written by Eric Young (eyay@cryptsoft.com).

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7) RZB / RCF obere Reihe

8) RZB / RCF untere Reihe

1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)

2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)

3) Observe the tube manufacturer's instructions.

4) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".

5) recommended for high-speed centrifugation

6) Load only each second position of the rotor

7) RZB / RCF top row

8) RZB / RCF bottom row